

Quick-Start Protocol

QIAseq® FastSelect™ –5S/16S/23S Kit with the TruSeq® Stranded Library Preparation

The QIAseq FastSelect –5S/16S/23S Kit (cat. nos. 335925, 335927, 335929) can be used for bacterial 5S/16S/23S rRNA removal with the TruSeq Stranded mRNA Library Prep (Illumina cat. no. 20020594, 20020595). Upon receipt, the FastSelect 5S/16S/23S tube should be immediately stored at –30 to –15°C in a constant-temperature freezer. All remaining components should immediately be stored in a refrigerator at 2–8°C.

Further information

- *QIAseq FastSelect –5S/16S/23S Handbook*: www.qiagen.com/HB-2695
- Safety Data Sheets: www.qiagen.com/safety
- Technical assistance: support.qiagen.com

Notes before starting

- The TruSeq Stranded mRNA Library Prep is required for use with this protocol.
Note: With this protocol, do not perform mRNA purification. Instead, follow the steps outlined below before proceeding to the designated step in the *TruSeq Stranded mRNA Reference Guide*. By doing this, a stranded total RNA library prep will be performed.
- For more information, please refer to the *TruSeq Stranded mRNA Reference Guide* (1000000040498).

Procedure

1. Thaw template RNA on ice. Gently mix, briefly centrifuge to collect residual liquid from the sides of the tubes, and then return to ice.



2. Prepare reagents required for RNA fragmentation and QIAseq FastSelect rRNA removal:
 - 2a. Bring FastSelect 5S/16S/23S, FastSelect FH Buffer, and Nuclease-free Water to room temperature.
 - 2b. **Important:** Only for FastSelect 5S/16S/23S, incubate the tube at 37°C for 5 min, and then vortex to dissolve the precipitate.
 - 2c. Mix all other reagents vigorously by vortexing, and then briefly centrifuge.
3. On ice, prepare the fragmentation/depletion reaction according to Table 1. Briefly centrifuge, mix by pipetting up and down 10 times, and then centrifuge briefly again.
4. Incubate as described in Table 2.

Important: Perform steps 2–9 regardless of the time and temperature in step 1.
5. Add 19.5 µl QIAseq Beads (prewarmed to room temperature and mixed by vortexing) to the 15 µl reaction. Mix thoroughly by vortexing, and incubate for 5 min at room temperature.
6. Centrifuge in a table-top centrifuge until the beads are completely pelleted (~2 min).
7. Place the tubes/plate on a magnetic rack for 2 min. Once the solution has cleared, with the beads still on the magnetic stand, carefully remove and discard the supernatant.

Important: Do not discard the beads, because they contain the RNA of interest.
8. Add 15 µl of Nuclease-free Water and 19.5 µl of QIAseq Bead Binding Buffer. Mix thoroughly by vortexing, and incubate for 5 min at room temperature.
9. Centrifuge in a table-top centrifuge until the beads are completely pelleted (~2 min).
10. Place the tubes/plate on a magnetic rack for 2 min. Once the solution has cleared, with the beads still on the magnetic stand, carefully remove and discard the supernatant.

Important: Do not discard the beads, because they contain the RNA of interest.
11. With the beads still on the magnetic stand, add 200 µl of 80% ethanol. Wait 30 s. Carefully remove and discard the wash.
12. Repeat the ethanol wash.

Important: Completely remove all traces of ethanol after this second wash. Remove the ethanol with a 200 µl pipette first, and then use a 10 µl pipette to remove any residual ethanol that will settle.

Table 1. Setup of fragmentation/depletion reactions

Component	Volume/reaction
Total RNA (100 ng – 1 µg)	Variable
FastSelect FH Buffer	1.5 µl
FastSelect 5S/16S/23S*	1 µl
Nuclease-free Water	Bring total reaction volume to 15 µl
Total volume	15 µl

* If needed, add 1 µl QIAseq FastSelect –rRNA HMR, 1 µl QIAseq FastSelect –Globin, 1 µl QIAseq FastSelect –rRNA Plant, and/or 1 µl QIAseq FastSelect –Yeast. To keep the total volume at 15 µl, correspondingly reduce the Nuclease-free Water by the volume of additional FastSelect products that have been added.

Table 2. Combined QIAseq fragmentation and FastSelect hybridization protocol

Note	Step	Mean insert size ~175–225 bp
Steps 2–9 are performed regardless of Input RNA quality. They need to be performed whether the RNA is high quality, moderate quality, FFPE, or degraded.	1*	8 min at 89°C*†
	2	2 min at 75°C
	3	2 min at 70°C
	4	2 min at 65°C
	5	2 min at 60°C
	6	2 min at 55°C
	7	2 min at 37°C
	8	2 min at 25°C
	9	Hold at 4°C

* For high-quality RNA (RIN ≥8), 8 min at 89°C is the basic recommendation. Alternatively, 4 min at 89°C provides an average insert size of ~325–375 bp, and 12 min at 89°C provides an average insert size of ~125–175 bp.

† If working with RNA with RIN values <8, please refer to *QIAseq FastSelect –5S/16S/23S Handbook*, Table 1, for fragmentation recommendations.

13. With the beads still on the magnetic stand, air-dry at room temperature for 6–10 min until all liquid has evaporated but without overdrying the beads.

Note: Visually inspect the pellet to confirm that it is completely dry.

14. Remove the beads from the magnetic stand, and elute the RNA from the beads by adding 19 µl PPF Buffer. Mix well by pipetting or vortexing, and then allow the tubes to sit at room temperature for 2 min to appropriately hydrate the beads.

Note: PPF Buffer is a component from the TruSeq Stranded mRNA Library Prep.

15. Return the tubes/plate to the magnetic rack until the solution has cleared.
16. Transfer 17 μ l of the supernatant to clean tubes/plate. Alternatively, the samples can be stored at -90 to -65°C in a constant-temperature freezer.
17. Refer to the *TruSeq Stranded mRNA Reference Guide* and immediately proceed to and perform "Synthesize First Strand cDNA".
Note: From the *TruSeq Stranded mRNA Reference Guide*, the procedural step "Place the RBP plate on the magnetic stand and wait until the liquid is clear (~5 min)" is not applicable.
18. Follow the *TruSeq Stranded mRNA Reference Guide* to perform all remaining library construction steps.
Important: It is highly recommended to dilute the Illumina adapters 2-fold compared to what is suggested in the default Illumina protocol.

Document Revision History

Date	Changes
07/2020	Clarified storage condition. Included 37°C incubation for FastSelect 5S/16S/23S as a default step in the procedure. Included options to include FastSelect $-$ rRNA Plant and/or FastSelect $-$ rRNA Yeast in the procedure. Replaced all precise storage temperatures with temperature ranges.



Scan QR code for handbook.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.

Trademarks: QIAGEN[®], Sample to Insight[®], QIAseq[®], FastSelect[™] (QIAGEN Group); TruSeq[®] (Illumina, Inc.). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

07/2020 HB-2706-003 © 2020 QIAGEN, all rights reserved.