

---

November 2015

# Investigator<sup>®</sup> STAR<sup>™</sup> Lyse&Prep Kit Handbook

For automated purification  
of DNA from forensic  
samples using STAR Q  
SP/AS instruments

---

# Contents

Kit Contents .....	3
Storage .....	4
Intended Use .....	4
Safety Information .....	5
Quality Control .....	5
Introduction .....	6
Principle and Procedure .....	6
Equipment and Reagents to Be Supplied by User .....	8
Important Notes .....	9
Starting material .....	9
Precipitate in Buffer ATL .....	9
Precipitate in Buffer QSL3 .....	9
Lysis with Proteinase K .....	10
Quantification of DNA .....	10
Protocol: Pretreatment for Various Casework and Reference Samples .....	12
Troubleshooting Guide .....	14
Ordering Information .....	15

# Kit Contents

<b>Investigator STAR Lyse&amp;Prep Kit</b>	<b>(400)</b>
<b>Catalog no.</b>	<b>931447</b>
<b>Number of preps</b>	<b>400</b>
Buffer ATL	5 x 50 ml
Buffer QSL3*	5 x 45 ml
Buffer QSW1*	5 x 50 ml
Buffer QSW2	5 x 160 ml
Bead Suspension G	5 x 15 ml
Buffer ATE	5 x 20 ml
Proteinase K	2 x 6 ml
Carrier RNA	2 x 310 µg
Q-Card†	1
Quick-Start Protocol	1
Handbook	1

\* Contains a guanidine salt. Not compatible with disinfectants containing bleach. See page 5 for Safety Information.

† The information encoded in the bar code on the Q-Card is needed for reagent data tracking.

---

## Storage

The Investigator STAR Lyse&Prep Kit is shipped at ambient temperature. All buffers and reagents can be stored at room temperature (15–25°C). Do not freeze the reagent bottles. When stored properly, the reagents are stable until the expiration date on the Q-Card. Lyophilized carrier RNA is stable until the expiration date on the Q-Card, when stored at room temperature.

The ready-to-use Proteinase K solution is stable for up to 1 year after delivery, when stored at room temperature.

## Intended Use

The Investigator STAR Lyse&Prep Kit is intended for molecular biology applications in forensic, human identity and paternity testing. This product is not intended for the diagnosis, prevention or treatment of a disease.

All due care and attention should be exercised in the handling of the products. We recommend all users of QIAGEN products to adhere to the NIH guidelines that have been developed for recombinant DNA experiments, or to other applicable guidelines.

# Safety Information

When working with chemicals, always wear a suitable lab coat, disposable gloves and protective goggles. For more information, please consult the appropriate safety data sheets (SDSs). These are available online in convenient and compact PDF format at [www.qiagen.com/safety](http://www.qiagen.com/safety) where you can find, view and print the SDS for each QIAGEN kit and kit component.



**CAUTION: DO NOT add bleach or acidic solutions directly to the sample preparation waste.**

Buffers QSL3 and QSW1 contain guanidine salts, which can form highly reactive compounds when combined with bleach. If liquid containing these buffers is spilt, clean with suitable laboratory detergent and water. If the spilt liquid contains potentially infectious agents, clean the affected area first with laboratory detergent and water, and then with 1% (v/v) sodium hypochlorite.

## Quality Control

In accordance with QIAGEN's ISO-certified Quality Management System, each lot of Investigator STAR Lyse&Prep Kits is tested against predetermined specifications to ensure consistent product quality.

Functional QC testing ensures that the Investigator STAR Lyse&Prep Kit meets the high standards required by forensic scientists.

---

# Introduction

The STAR Q SP/AS instrument and the Investigator STAR Lyse&Prep Kit reproducibly automate purification of genomic DNA, encountered in forensic and human identity applications. Purification is efficient, and purified DNA performs well in downstream analyses, such as quantitative PCR and STR analysis.

## Principle and Procedure

Magnetic-particle technology combines the speed and efficiency of silica-based DNA purification with the convenient handling of magnetic particles. DNA is isolated from lysates in one step through its binding to the silica surface of the particles in the presence of a chaotropic salt. The particles are separated from the lysates using a magnet. The DNA is then efficiently washed and eluted in modified TE buffer (Buffer ATE). The user can choose elution volumes of 50 or 100  $\mu$ l.

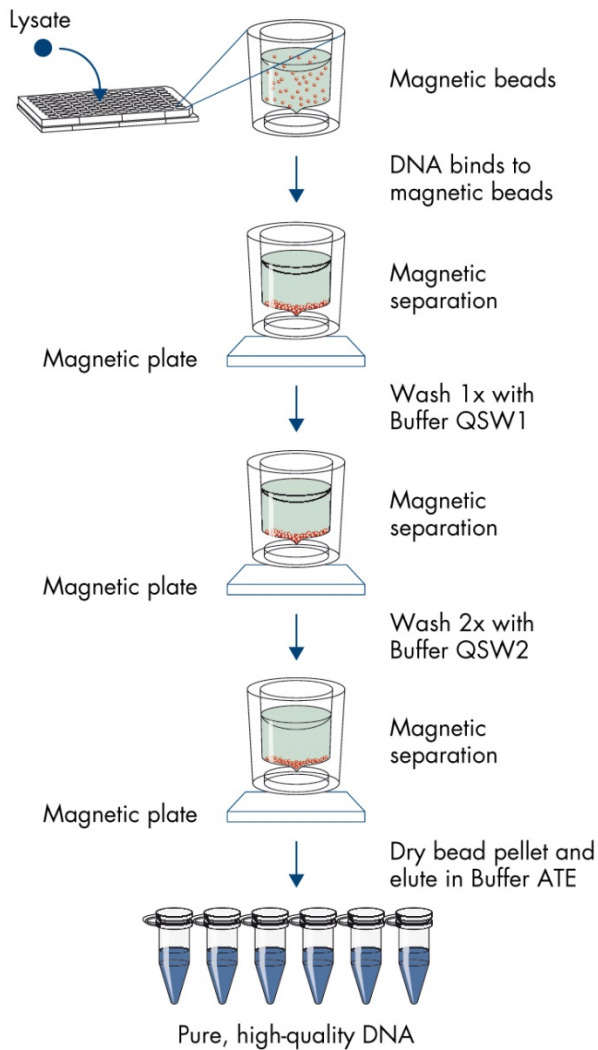


Figure 1. The STAR Q SP/AS workflow

# Equipment and Reagents to Be Supplied by User

When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, consult the appropriate safety data sheets (SDSs), available from the product supplier.

## All protocols

- Sample and elution labware according to Table 1
- Thermomixer
- Centrifuge
- Vortexer
- Pipets and pipet tips (to prevent cross-contamination, we strongly recommend the use of pipet tips with aerosol barriers)

For purification of DNA from semen stains or hair

- 1 M dithiothreitol (DTT)

**Table 1. Labware for sample input and elution**

Manufacturer	Type	Catalogue no.	Use
Treff*	2 ml	96.09329.9.01	Sample input/elution
Eppendorf®	1.5 ml Safe-Lock	0017 010.417-03/1109	Sample input/elution
Eppendorf	2 ml Safe-Lock	0017 010.425-03/1109	Sample input/elution
Sarstedt®	1.5 ml Safe-Lock	72.706	Sample input/elution
Sarstedt	2 ml Safe-Lock	72.695.500	Sample input/elution
Sarstedt	1.5 ml	72.690.001	Sample input/elution
Sarstedt	2 ml	72.691	Sample input/elution
VWR®	96-Well-Plate, 2.2 ml	732-0585	Sample input

\* Collection tube of the Investigator Lyse&Spin Basket Kit.



---

# Important Notes

## Starting material

The amount of starting material for use in Investigator STAR Lyse&Prep Kit procedures can vary greatly, depending on the amount of DNA in the sample. Specific guidance for starting amounts is given in the individual protocols, and Table 2.

## Purification of low amounts of DNA

For purification of DNA from very small amounts of sample, such as touch DNA samples, we recommend adding carrier RNA. For samples containing larger amounts of DNA, addition of carrier RNA is optional. Add 310  $\mu\text{l}$  TE buffer or water to the tube containing 310  $\mu\text{g}$  lyophilized carrier RNA to obtain a solution of 1  $\mu\text{g}/\mu\text{l}$ . Dissolve the carrier RNA thoroughly, divide in to conveniently sized aliquots, and store at  $-20^{\circ}\text{C}$ . Do not freeze-thaw the aliquot of carrier RNA more than 3 times. Carrier RNA should be added to the sample after the lysis is completed to avoid degradation. Carrier RNA can be added to the binding buffer prior to setting up the instrument run. Use the mixture of binding buffer and carrier RNA within the same day.

## Precipitate in Buffer ATL

Before starting the procedure, check whether precipitate has formed in Buffer ATL. If necessary, dissolve by heating to  $70^{\circ}\text{C}$  with gentle agitation.

## Precipitate in Buffer QSL3

Buffer QSL3 may form a precipitate upon storage. If necessary, redissolve by mild agitation at  $37^{\circ}\text{C}$  and then place at room temperature ( $15$ – $25^{\circ}\text{C}$ ).

---

## Lysis with Proteinase K

The Investigator STAR Lyse&Prep Kit contains Proteinase K, which is the enzyme of choice for lysis buffers used in Investigator STAR Lyse&Prep protocols. Proteinase K is a recombinant protein expressed in *Pichia pastoris*, and is particularly suitable for short digestion times. It possesses a high specific activity and remains stable over a wide range of temperatures and pH values, with substantially increased activity at higher temperatures. The activity of the Proteinase K solution is 600 mAU/ml solution (or 40 mAU/mg protein).

## Quantification of DNA

Degraded, inhibited or mixed DNA samples are common in forensic casework and other human identity testing applications. Such samples can create challenges in STR analysis. Prior quantification of the purified DNA using real-time PCR is recommended and reduces the need to repeat downstream analyses. This greatly reduces costs and time, and improves the statistical relevance of results. Investigator Quantiplex® Kits, use quantitative real-time PCR to quantify human DNA in a sample. These kits also detect if a sample contains sufficient DNA to enable DNA fingerprinting analysis and if the sample contains inhibitors that may interfere with downstream applications. They are designed for use in a whole range of human identity and forensics applications, and are available in 2 versions: 1) the Investigator Quantiplex Kit, for quantification of total human DNA, and 2) the Investigator Quantiplex HYres Kit, for quantification of total human and human male DNA.

**Table 2. Protocol information for different sample types**

<b>Sample type</b>	<b>Sample amount</b>	<b>Buffer ATL</b>	<b>Proteinase K</b>	<b>DTT (1 M)</b>
Blood/saliva	Up to 50 µl	280 µl/475 µl*	20 µl/25 µl*	–
Surface swabs	1 swab	280 µl/475 µl*	20 µl/25 µl*	–
Chewing gum	Up to 40 mg	280 µl/475 µl*	20 µl/25 µl*	–
Cigarette butts	1 cm <sup>2</sup>	280 µl/475 µl*	20 µl/25 µl*	–
Paper/similar materials	0.5–2.5 cm <sup>2</sup>	280 µl/475 µl*	20 µl/25 µl*	–
Nail scrapings	Up to 40 mg	260 µl/455 µl*	20 µl/25 µl*	20 µl
Nail clippings	1	260 µl/455 µl*	20 µl/25 µl*	20 µl
Hair	0.5–1 cm	260 µl/455 µl*	20 µl/25 µl*	20 µl
Tissues	Up to 10 mg	280 µl/475 µl*	20 µl/25 µl*	–
Blood or saliva stains	0.5 cm <sup>2</sup>	280 µl/475 µl*	20 µl/25 µl*	–
Semen stains	0.5 cm <sup>2</sup>	260 µl/455 µl*	20 µl/25 µl*	20 µl
Buccal swabs	1 swab	280 µl/475 µl*	20 µl/25 µl*	–

\* If using the Investigator Lyse&Spin Basket Kit.

---

# Protocol: Pretreatment for Various Casework and Reference Samples

This protocol is designed for isolation of total DNA (genomic and mitochondrial) from various types of casework and reference samples. The protocol describes the preliminary lysis using Proteinase K.

## Notes before starting

- Before beginning the procedure, read “Important Notes” on page 9.
- We recommend using the Investigator Lyse&Spin Basket Kit when solid sample materials have to be removed from the lysate. If using this kit, please follow the Pretreatment protocol, as outlined under “Procedure using the Investigator Lyse&Spin Basket Kit” (page 13). The Lyse&Spin Basket Kit collection tubes can be used as sample tubes.
- Heat a thermomixer to 56°C for the Proteinase K digest in step 3.

## Procedure using 300 µl lysis volume

1. Place the sample into a sample tube or deep-well block.
2. Set up the Proteinase K digest according to the information provided in Table 2, page 11. Mix sample thoroughly by vortexing for 10 s.
3. Incubate at 56°C for 15 min to overnight in a thermomixer, shaking at 900 rpm.  
Incubation for 15 min may be sufficient to recover adequate DNA for STR typing from samples containing abundant DNA. More than 1 h is recommended where a low amount of DNA is expected. When using in a deep-well block, ensure proper sealing with an adhesive tape.
4. After incubation, perform a brief centrifugation to collect liquid at the bottom of the tubes or the S-Block.
5. Transfer 300 µl lysate to a new tube or deep-well block.

---

**Take care not to transfer any solid sample materials. See Table 1 for suitable labware.**

6. Continue with DNA Purification on the STAR Q SP/AS instrument.

**The software guides you through the worktable setup. Choose 300  $\mu$ l as lysate volume.**

#### Procedure using the Investigator Lyse&Spin Basket Kit

1. Place the stained material into the Investigator Lyse&Spin Basket, inserted into a 2 ml microcentrifuge tube (provided).
2. Set up the Proteinase K digest according to the information provided in Table 2. Mix sample thoroughly by vortexing for 10 s.

**Note:** The Investigator Lyse&Spin Basket requires 500  $\mu$ l lysate volume.

3. Incubate at 56°C for 1 h to overnight in a thermomixer, shaking at 900 rpm.
4. Centrifuge for 1 min at a minimum of 10,000  $\times$  g.

**Note:** Keep the lid closed during centrifugation.

**Note:** Up to 20,000  $\times$  g can be used for centrifugation.

**Note:** Make sure that no liquid remains in the basket after centrifugation. If necessary, repeat the centrifugation until all liquid has passed through the membrane. If larger pieces of chewing gum are processed, clogging of the basket can be avoided by pressing the chewing gum against the sides of the basket.

5. Discard the basket including the solid sample substrate.
6. Continue with DNA Purification on the STAR Q SP/AS instrument.

**The software guides you through the worktable setup. Choose 500  $\mu$ l as lysate volume.**

# Troubleshooting Guide

This troubleshooting guide may be helpful in solving any problems that may arise. For more information, see also the Frequently Asked Questions page at our Technical Support Center: [www.qiagen.com/FAQ/FAQList.aspx](http://www.qiagen.com/FAQ/FAQList.aspx). The scientists in QIAGEN Technical Services are always happy to answer any questions you may have about either the information and/or protocols in this handbook or sample and assay technologies (for contact information, visit [www.qiagen.com](http://www.qiagen.com)).

## Comments and suggestions

---

### Low DNA yield

- |    |                                       |   |
|----|---------------------------------------|---|
| a) | Insufficient lysis                    | Increase lysis time.  |
| b) | Insufficient sample aspirated         | Check sample input for remaining lysate.  |
| c) | Insufficient amount of magnetic beads | Carefully mix bead suspension before transferring to tubes to load on the instrument. |

### DNA does not perform well in downstream applications

- |    |  |   |
|----|--|---|
| a) | Insufficient DNA used in downstream applications | If possible, repeat the downstream application using more eluate.   |
| b) | Excess of magnetic beads                         | Excess of magnetic beads may lead to insufficient drying before elution, resulting in ethanol carry over. Carefully mix bead suspension before transferring to tubes to load on the instrument. |

# Ordering Information

<b>Product</b>	<b>Contents</b>	<b>Cat. no.</b>
Investigator STAR Lyse&Prep Kit (400)	For 400 preps of 300 µl each from casework and reference samples	931447
STAR Q SP/AS Instrument	Liquid-handling instrument, PC, STAR Q SP/AS software	9002650
STAR Q Lyse&Prep Instrument	Liquid-handling instrument, PC, STAR Q Lyse&Prep software	9002653
Investigator Lyse&Spin Basket Kit (50)	50 pouches containing 50 baskets and 100 collection tubes	19597
Investigator Lyse&Spin Basket Kit (250)	10 pouches containing 5 x 50 baskets and 5 x 50 collection tubes	19598

---

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at [www.qiagen.com](http://www.qiagen.com) or can be requested from QIAGEN Technical Services or your local distributor.

Visit [www.qiagen.com](http://www.qiagen.com) to find out more!



Trademarks: QIAGEN<sup>®</sup>, Sample to Insight<sup>®</sup>, Investigator<sup>®</sup>, Quantiplex<sup>®</sup> (QIAGEN Group); Eppendorf<sup>®</sup> (Eppendorf AG); STAR<sup>™</sup> (Hamilton Company); Sarstedt<sup>®</sup> (Sarstedt AG and Co.); VWR<sup>®</sup> (VWR International, LLC). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

#### **Limited License Agreement for Investigator STAR Lyse&Prep Kit**

Use of this product signifies the agreement of any purchaser or user of the product to the following terms:

1. The product may be used solely in accordance with the protocols provided with the product and this handbook and for use with components contained in the kit only. QIAGEN grants no license under any of its intellectual property to use or incorporate the enclosed components of this kit with any components not included within this kit except as described in the protocols provided with the product, this handbook, and additional protocols available at [www.qiagen.com](http://www.qiagen.com). Some of these additional protocols have been provided by QIAGEN users for QIAGEN users. These protocols have not been thoroughly tested or optimized by QIAGEN. QIAGEN neither guarantees them nor warrants that they do not infringe the rights of third-parties.
2. Other than expressly stated licenses, QIAGEN makes no warranty that this kit and/or its use(s) do not infringe the rights of third-parties.
3. This kit and its components are licensed for one-time use and may not be reused, refurbished, or resold.
4. QIAGEN specifically disclaims any other licenses, expressed or implied other than those expressly stated.
5. The purchaser and user of the kit agree not to take or permit anyone else to take any steps that could lead to or facilitate any acts prohibited above. QIAGEN may enforce the prohibitions of this Limited License Agreement in any Court, and shall recover all its investigative and Court costs, including attorney fees, in any action to enforce this Limited License Agreement or any of its intellectual property rights relating to the kit and/or its components.

For updated license terms, see [www.qiagen.com](http://www.qiagen.com).

1099541 HB-1983-002 11/2015 © 2015

---

## Notes

---

## Notes

---

Ordering [www.qiagen.com/contact](http://www.qiagen.com/contact) | Technical Support [support.qiagen.com](http://support.qiagen.com) | Website [www.qiagen.com](http://www.qiagen.com)