

QIAseq® 16S/ITS Screening Panel

Further information

- *QIAseq 16S/ITS Panels Handbook*: www.qiagen.com/HB-2547
- Safety Data Sheets: www.qiagen.com/safety
- Technical assistance: toll-free 00800-22-44-6000, or www.qiagen.com/contact

Important points before starting

- Use high quality DNA.
- If DNA concentration is >1 ng/ μ l, then dilute to 1 ng/ μ l. If DNA concentration is <1 ng/ μ l, then use 1 μ l of sample per PCR reaction.
- Ensure reactions are thoroughly mixed, as well as prepared and incubated at the recommended temperatures.
- 80% Ethanol should be freshly prepared using nuclease-free water and mixed thoroughly by vortexing.
- QIAseq Beads need to be homogenous. Thoroughly resuspend the beads immediately before use and process the beads quickly. If a delay in the protocol occurs, simply vortex the beads again.

Preparation of QIAseq 16S/ITS Screening Panel PCR Reaction

1. On ice, prepare three PCR reactions per gDNA sample by following Table 1. Mix well by pipetting up and down 10 times or by vortexing, then spin down.

Table 1. Preparation of QIAseq 16S/ITS Region Panel PCR

Component	Per Sample		
	Panel Pool 1	Panel Pool 2	Panel Pool 3
Microbial DNA Sample	1 µl	1 µl	1 µl
UCP Multiplex Master Mix	2.5 µl	2.5 µl	2.5 µl
Panel Pool 1	1 µl	—	—
Panel Pool 2	—	1 µl	—
Panel Pool 3	—	—	1 µl
UCP PCR Water	5.5 µl	5.5 µl	5.5 µl
Total Volume	10 µl	10 µl	10 µl

2. Incubate the reactions in a thermal cycler as described in Table 2.

Table 2. Setup of QIAseq 16S/ITS PCR Reaction

Step	Time	Temperature
Hold	2 mins	95°C
3-step cycling		
Denaturation	30 secs	95°C
Annealing	30 secs	50°C
Extension	2 mins	72°C
12 cycles*		
Final Extension (1 cycle)	7 mins	72°C
Hold	∞	4°C

*If samples are low biomass or contain low abundance of bacteria, use 20 cycles of 16S PCR.

3. After PCR is finished, add 20 µl of UCP Water to each of the PCR reactions.
4. Pool the PCR reactions into a single LoBind tube or well of a 96-well PCR plate.

-
5. Add 99 μ l of QIAseq Beads and mix well by pipetting up and down 12 times.
 6. Incubate for 5 minutes at room temperature.
 7. Place the tubes/plate on a magnetic rack for 5 minutes and discard the supernatant.
 8. Centrifuge and carefully remove any residual liquid. Add 105 μ l of nuclease-free water. Mix by pipetting up and down 12 times until the beads are fully resuspended.
 9. Place the tubes/plate on a magnetic rack for 5 minutes. Carefully transfer 100 μ l of the supernatant that contains the 16S/ITS PCR product to new tubes or plate.
 10. Add 110 μ l of QIAseq Beads to each sample. Mix well by pipetting up and down.
 11. Incubate for 5 minutes at room temperature.
 12. Place the tubes/plate on a magnetic rack for 5 minutes. After the solution has cleared, carefully remove and discard the supernatant.
 13. Wash with 200 μ l of 80% ethanol. Carefully remove and discard the wash.
 14. Repeat the ethanol wash and ensure that all traces of ethanol have been removed.
 15. With the tubes/plate still on the magnet, air dry at room temperature for 10 minutes.
 16. Elute the DNA from the beads by adding 35 μ l UCP PCR Water. Mix well.
 17. Return the tubes/plate to the magnetic rack until the solution has cleared.
 18. Transfer 32.5 μ l of the supernatant to clean tubes/plate.
 19. Proceed to Protocol: *Preparation of QIAseq 16S/ITS Screening Panel Sample Index PCR Reaction*.
-

Preparation of QIAseq 16S/ITS Screening Panel Sample Index PCR Reaction

1. To the tubes/plate containing the 16S/ITS PCR Product, add the components according to Table 3. Mix well by pipetting up and down 12 times or by vortexing, then spin down.

Table 3. Preparation of QIAseq 16S/ITS Sample Index PCR Reaction

Step	Per Sample	HT Array
16S/ITS Region Panel PCR Product	32.5 µl	32.5 µl
UCP Master Mix	12.5 µl	12.5 µl
p5-RS2-ID# (4 µM) *	2.5 µl	-
p7-FS2-ID# (4 µM) *	2.5 µl	-
UCP Water		5 µl
Total Volume	50 µl	50 µl

*Use a unique p5-RS2-ID# + p7-FS2-ID# combination for each Microbial DNA Sample.

2. Incubate the reactions in a thermal cycler according to Table 4.

Table 4. Setup of QIAseq 16S/ITS Sample Index PCR Reaction

Step	Time	Temperature
Hold	2 mins	95°C
3-step cycling		
Denaturation	30 secs	95°C
Annealing	30 secs	60°C
Extension	2 mins	72°C
14 cycles		
Final Extension (1 cycle)	7 mins	72°C
Hold	∞	4°C

3. Remove the tubes/plate from the thermal cycler and briefly centrifuge.

-
4. Add 45 µl of QIAseq Beads and mix well by pipetting up and down 12 times.
 5. Incubate for 5 min at room temperature.
 6. Place the tubes/plate on a magnetic rack for 5 minutes and discard the supernatant.
 7. Wash with 200 µl of 80% ethanol. Carefully remove and discard the wash.
 8. Repeat the ethanol wash and ensure that all traces of ethanol have been removed.
 9. With the tubes/plate still on the magnet, air dry at room temperature for 10 minutes.
 10. Elute the DNA from the beads by adding 30 µl nuclease-free water. Mix well.
 11. Return the tubes/plate to the magnetic rack until the solution has cleared.
 12. Transfer 25 µl of the supernatant to clean tubes/plate.
 13. Proceed to *Protocol: Library QC and Quantification*. Alternatively, the completed QIAseq 16S/ITS Screening Panel Sequencing Library can be stored at -20°C in a constant-temperature freezer.
-



Scan QR code for handbook.

Trademarks: QIAGEN®, Sample to Insight®, QIAseq® (QIAGEN Group); LoBind® (Eppendorf AG); Ion Torrent®, (Thermo Fisher Scientific or its subsidiaries).
For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.
1115002 12/2018 HB-2565-002

© 2018 QIAGEN, all rights reserved.

Ordering www.qiagen.com/shop | Technical Support support.qiagen.com | Website www.qiagen.com