

Protocol Sheet

Applied Biosystems® 7900HT real-time PCR run setup instructions for miScript® miRNA PCR Arrays

Important points before starting

- Please read the handbook supplied with the miScript miRNA PCR Array, paying careful attention to the “Safety Information” and “Important Notes” sections, before beginning this procedure.
- Please make sure the real-time PCR instrument is working properly. Refer to the manufacturer’s Installation and Maintenance manual if needed.

Procedure

Creation of PCR protocol template

1. Open the ABI 7900HT SDS Version 2.4 software on the desktop of the computer that is connected to the ABI 7900HT instrument.
2. Select “File” > “New.” The “New Document Wizard” dialog box will appear.
 - Select “Assay”: “Standard Curve (AQ).”
 - Select “Container”: “384-well Clear Plate.”
 - Select “Template”: “Blank Template.”
 - The field for “Barcode” is optional.
 - Click “OK.”
 - Under the “Setup” tab, click the “Add Detector...” button at the bottom of the tab.
 - In the “Detector Manager” window, select SYBR® for the reporter dye, click the “Copy To Plate Document” button, and click the “Done” button (see Figure 1). Ensure that ROX is selected for “Passive Reference.”



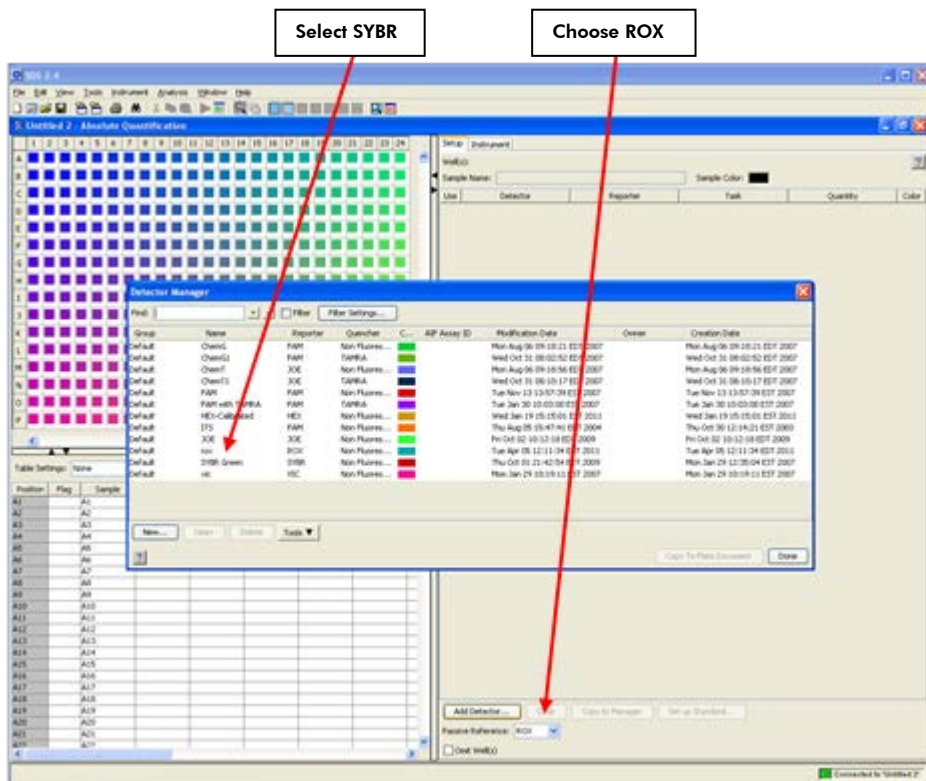


Figure 1. "Detector Manager" screen.

- Click the square button in the upper left corner of the 384-well layout panel (between the letter 'A' and '1') to select all wells. Once selected, the 384-well table will be highlighted in yellow. Click the box next to SYBR Green. This selects SYBR Green as the detector for all wells.

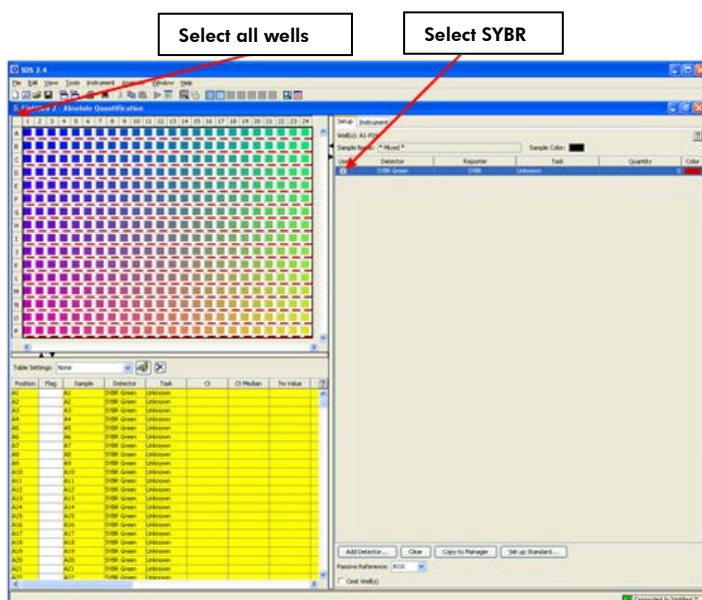


Figure 2. "Absolute Quantification" screen.

- Click the “Instrument” tab, then the “Thermal Cycler” > “Thermal Profile” tab (see Figure 3).
 - Highlight and delete Stage 1 (50°C for 2:00 minutes) by clicking the mouse, dragging it across Stage 1, and clicking the “Delete Step” button.
 - Stage 1: Enter 95.0°C for 15:00 minutes.
 - Stage 2 (3 Steps)
 - Step 1: 94.0°C for 0:15 (15 seconds)
 - Step 2: 55.0°C for 0:30 (30 seconds)
 - Step 3: 70.0°C for 0:30 (30 seconds)
 - Enter 40 for “Repeats.”
 - Click the “Add Dissociation Stage” button. This will add the pre-set dissociation stage as Stage 3.
 - Ensure that “Standard” is checked for “Mode.”
 - Type 10 for “Sample Volume (µl).”

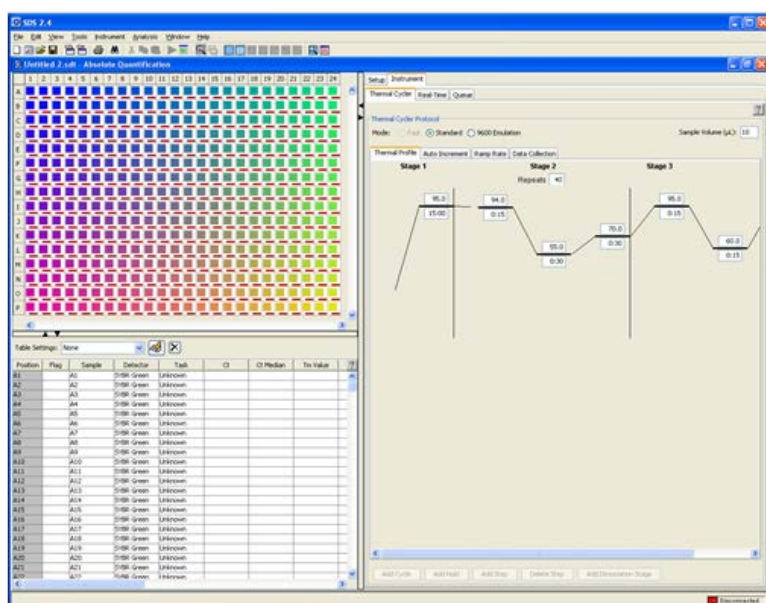


Figure 3. Thermal profile.

- Select “File” > “Save As” to save the template file. Save the file as SDS Templates (*.sdt) with the filename “miScript_miRNA_PCR_Array_Template” (click “Save”).

Note: Alternatively, download the PCR protocol template file

(miScript_miRNA_Template_ABI7900_M) from the QIAGEN website, www.qiagen.com.

Performing real-time PCR detection

3. If the thermocycler is off, press the power button to switch on the instrument. Wait for the instrument to boot and display the "Power" status light. Switch on the computer connected to the thermocycler.
4. Ensure that the plate has been centrifuged for 1 min. at 1000 g to remove any bubbles.
5. Open the ABI 7900HT SDS Version 2.4 software.
6. Select "File" > "New." In the "New Document Wizard" dialog box, select "Browse" to load the "miScript_miRNA_PCR_Array_Template" file. Then click "Finish." This will load the previously saved setup to the new plate document.
7. Save the new document under a new filename as an SDS document (*.sds).
8. Click the "Instrument" tab. Click "Real-Time" tab, then click "Connect to Instrument" to connect the computer to the thermal cycler. Click "Open/Close" to open the plate tray and place your plate in the precision plate holder with A1 in the top left corner. Click "Start Run" to begin the PCR run. Wait 30 seconds to 1 min. for the initial priming, and then the run should start. The estimated run time will then appear on the screen.

After the PCR run

9. When the PCR run is complete, a small dialog box stating "The run completed successfully" will appear on the screen. Click "OK"; this will close the box.
10. To determine C_T values, set Baseline and Threshold as described in the *miScript miRNA PCR Array Handbook*.
11. The C_T values will be displayed in the lower left panel for each well. To export the result to an Excel® spreadsheet, select "File" > "Export" > "Results Table." The file will be saved as a tab-delimited text file, and can be opened using Microsoft® Excel.

The miScript miRNA PCR Array is intended for molecular biology applications. This product is not intended for the diagnosis, prevention, or treatment of a disease.

For up-to-date licensing and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN handbooks can be requested from QIAGEN Technical Service or your local QIAGEN distributor. Selected handbooks can be downloaded from www.qiagen.com/literature. Safety data sheets (SDS) for any QIAGEN product can be downloaded from www.qiagen.com/safety.

Trademarks: miScript®, QIAGEN® (QIAGEN Group); Applied Biosystems®, SYBR® (Life Technologies Corporation); Microsoft®, Excel® (Microsoft Corporation).
Dec-12 © 2012 QIAGEN, all rights reserved.

www.qiagen.com

Australia ■ 1-800-243-800

Austria ■ 0800/281010

Belgium ■ 0800-79612

Canada ■ 800-572-9613

China ■ 021-51345678

Denmark ■ 80-885945

Finland ■ 0800-914416

France ■ 01-60-920-930

Germany ■ 02103-29-12000

Hong Kong ■ 800 933 965

Ireland ■ 1800 555 049

Italy ■ 800 787980

Japan ■ 03-6890-7300

Korea (South) ■ 1544 7145

Luxembourg ■ 8002 2076

The Netherlands ■ 0800 0229592

Norway ■ 800-18859

Singapore ■ 65-67775366

Spain ■ 91-630-7050

Sweden ■ 020-790282

Switzerland ■ 055-254-22-11

UK ■ 01293-422-911

USA ■ 800-426-8157



Sample & Assay Technologies