

August 2023

Release Note: EZ2[®] Connect MDx Protocol Package v8

Dear valued customer,

At QIAGEN[®], we are continuously making improvements to provide the highest level of performance with our instruments and protocols. Consequently, we are pleased to inform you about the opportunity to upgrade your EZ2 Connect MDx instrument to Protocol Package v8. The protocol package update has no influence on the IVD protocols of EZ2 Connect MDx. If you are using protocols below, we highly recommend the protocol package update to further reduce the risks and to benefit from the improvements as detailed below.

Protocol name	Cat. no. of the kit	New protocol version	Changes to protocol package v7
EZ1&2 [™] DNA Blood 200 µL Kit	951034	4	More accurate run time. Additional measures to reduce the risk of lifting the tube up.
EZ1&2 DNA Blood 350 µL Kit	951054	5	More accurate run time. Sample volume displayed in the loading instructions.
EZ2 AdnaTest CTC Select Kit	395692	4	Additional measures to reduce the risk of heater time out with unstable or high environmental temperature.
EZ1&2 DNA FFPE Kit	954404, 954414	3	Additional measures to reduce the risk of splashing and mixing issues.
EZ2 RNA FFPE Kit	959734	4	Additional measures to reduce the risk of splashing and mixing issues. Sample volume displayed in the loading instructions.
EZ2 AllPrep [®] DNA/RNA FFPE Kit	954734	4	Additional measures to reduce the risk of splashing and mixing issues. More flexible choice of elution volumes.
EZ1&2 ccfDNA Kit	9027011	5	Additional measures to reduce the risk of heater time out with unstable or high environmental temperature. More flexible choice of elution volumes. New protocol for urine as starting material.



If you have further questions, please contact your local QIAGEN representative or contact our Technical Support Center at www.qiagen.com/support/technical-support.

Best regards,

Your EZ2 Connect Team