

Important Note

Use of Abnova™ RNA Marker Low with QIAxcel® ScreenGel®

Dear valued QIAxcel Connect customer,

For QIAxcel ScreenGel 2.1 and later, we are offering a Size Marker file to analyze RNA fragments of 20–500 nucleotides (available for download from www.qiagen.com/qiaxcel-connect-system). The Abnova RNA Marker Low (cat. no. R0001) is to be used in a concentration of 3.5 ng/μL (1:200 dilution with QX RNA HS Dilution Buffer), in combination with the QX RNA High Sensitivity Cartridge, QX RNA HS Alignment Marker, and ScreenGel Method Small RNA High-Sensitivity (available as web download from the QIAxcel webpage).

If you have further questions, please contact QIAGEN Technical Support. Visit support.qiagen.com for additional resources and contact information.

Best regards,

QIAGEN

Trademarks: QIAGEN®, Sample to Insight®, QIAxcel®, ScreenGel® (QIAGEN Group). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

06/2026 HB-3853-001 © 2026 QIAGEN, all rights reserved.