

# QIAGEN Supplementary Protocol

---

## Purification of REPLI-g<sup>®</sup> amplified DNA by LiCl/EtOH precipitation

This protocol is designed for the purification of 5–40 µg DNA amplified using the REPLI-g Single Cell Kit (cat. nos. 150343 and 150345), REPLI-g WTA Single Cell Kit (cat. nos. 150063 and 150065), or the REPLI-g Cell WGA & WTA Kit (cat. nos. 150052 and 150054) by lithium chloride/ethanol precipitation.

### Product use limitations

REPLI-g Kits are intended for molecular biology applications. These products are not intended for the diagnosis, prevention, or treatment of a disease.

All due care and attention should be exercised in the handling of the products. We recommend all users of QIAGEN<sup>®</sup> products to adhere to the NIH guidelines that have been developed for recombinant DNA experiments, or to other applicable guidelines.

### Equipment and reagents to be supplied by user

When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, consult the appropriate safety data sheets (SDSs), available from the product supplier.

- 1.5 or 2 ml microcentrifuge tubes
- Microcentrifuge
- Pipet tips (pipet tips with aerosol barriers for preventing cross-contamination are recommended)
- TE buffer (10 mM Tris-Cl; 1 mM EDTA, pH 8.0)
- Ethanol (96–100%)
- Ethanol (70%)
- Lithium chloride (7.5 M)
- EDTA (0.5 M, pH 8.0)



## Important points before starting

- All centrifugation steps should be performed at room temperature (15–25°C).
- Wear gloves throughout the entire procedure. In case of contact between gloves and sample, change gloves immediately.

## Procedure

- 1. After the REPLI-g reaction, equilibrate REPLI-g amplified DNA to room temperature (15–25°C) for no longer than 10 min.**

**Note:** If not purifying the complete volume, transfer the amplified DNA to a new 1.5 ml microcentrifuge tube.
- 2. Adjust the volume to 60 µl by adding the appropriate volume of TE buffer if necessary.**
- 3. Add 30 µl 7.5 M lithium chloride and 6 µl EDTA (0.5 M; pH 8.0) to REPLI-g amplified DNA, mix the sample by flicking the tube several times, centrifuge briefly, and let the sample incubate for 30 min at room temperature (15–25°C).**
- 4. Add 150 µl ethanol (96–100%). Mix carefully by flicking the tube several times.**

**Note:** Do not mix by pipetting up and down.

**Note:** A cotton-like precipitate will be visible.
- 5. Centrifuge at maximum speed (15000 x g) for 2 min. With most centrifuges, 15000 x g corresponds to 13000 rpm.**

**Note:** Do not centrifuge for more than 2 min. Centrifugation for longer periods will condense the precipitate, making it more difficult to dissolve.
- 6. Aspirate supernatant carefully using a pipet and discard the supernatant.**
- 7. Add 100 µl ethanol (70%) to the pellet. Flick the tube several times.**
- 8. Centrifuge at maximum speed for 1 min.**

**Note:** Do not centrifuge for more than 1 min. Centrifugation for longer periods will condense the precipitate, making it more difficult to dissolve.
- 9. Repeat steps 6–8.**
- 10. Aspirate supernatant carefully using a pipet and discard the supernatant.**
- 11. Centrifuge briefly to collect residual supernatant at the bottom of the tube.**
- 12. Aspirate residual supernatant carefully using a pipet and discard the supernatant.**

**Note:** Avoid aspiration of the pellet.
- 13. Incubate the microcentrifuge tube containing the precipitate of REPLI-g amplified DNA for 10 min at room temperature (15–25°C).**
- 14. Add 60 µl 1x TE buffer (pH 8.0) to dissolve REPLI-g amplified DNA.**

**Note:** If the DNA is not dissolved completely, dissolve the pellet overnight at 4–10°C.

**Note:** Mix carefully to avoid shearing of REPLI-g amplified DNA.

Purified DNA can be directly used for all downstream applications. It does not contain tags or labeled molecules. Concentration of purified DNA/cDNA is typically above 150 ng/μl if DNA/cDNA of a complete reaction is purified.

- 15. If not being used directly, store the amplified DNA/cDNA at –15 to –30°C until required for downstream applications. We recommend storage of the amplified DNA/cDNA at a concentration of at least 100 ng/μl.**

Amplified DNA/cDNA behaves like purified genomic DNA and has an approximate length of 2000 bp up to 70,000 bp. It is highly suited for use in a variety of downstream applications, particularly next-generation sequencing and quantitative PCR. See Table 1 for information on handling of amplified DNA/cDNA.

Optical density (OD) measurements can accurately determine the concentration of purified REPLI-g amplified DNA, for example, using the QIAxpert (QIAGEN, cat. no. 9002340). For applications where accurate quantification of double-strand DNA is especially important, such as using the Nextera library prep protocol to prepare an NGS library, concentration determination has to be performed using Quant-iT™ PicoGreen® dsDNA reagent (Life Technologies, cat no. P7581) or the Qubit® dsDNA BR Assay system (Life Technologies, cat. no. Q32850).

**Table 1. Applications and handling**

Downstream application	Use of amplified DNA/cDNA	QIAGEN products
NGS	Covaris®: 500 – 2000 ng* Nextera Library prep kit: 50 ng† GeneRead™ DNA Library Prep Kits†: 50–1000 ng fragmented DNA	GeneRead DNA Library Prep Kit
Real-time PCR, PCR	10 ng	QuantiTect®, QuantiFast®, QuantiNova™ Kits
Microarray	See supplier's instructions	–
Sanger sequencing, Pyrosequencing®	PCR has to be performed from the region of interest prior to sequencing. See advice for PCR.	PyroMark® products

\* Dependent on sequencing platform used.

† See dedicated supplementary protocols on [www.qiagen.com](http://www.qiagen.com).

QIAGEN handbooks can be requested from QIAGEN Technical Service or your local QIAGEN distributor. Selected handbooks can be downloaded from [www.qiagen.com/literature](http://www.qiagen.com/literature).

Safety data sheets (SDS) for any QIAGEN product can be downloaded from [www.qiagen.com/safety](http://www.qiagen.com/safety).

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at [www.qiagen.com](http://www.qiagen.com) or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, GeneRead™, PyroMark®, Pyrosequencing®, QuantiFast®, QuantiNova™, QuantiTect®, REPLI-g® (QIAGEN Group); Covaris® (Covaris, Inc.); PicoGreen®, QuantiT™ Qubit® (Life Technologies Corporation).

RG33 June-14 © 2014 QIAGEN, all rights reserved.

**www.qiagen.com**

**Australia** ■ 1-800-243-800

**Austria** ■ 0800/281010

**Belgium** ■ 0800-79612

**Canada** ■ 800-572-9613

**China** ■ 021-51345678

**Denmark** ■ 80-885945

**Finland** ■ 0800-914416

**France** ■ 01-60-920-930

**Germany** ■ 02103-29-12000

**Hong Kong** ■ 800 933 965

**Ireland** ■ 1800 555 049

**Italy** ■ 800 787980

**Japan** ■ 03-5547-0811

**Korea (South)** ■ 1544 7145

**Luxembourg** ■ 8002 2076

**The Netherlands** ■ 0800 0229592

**Norway** ■ 800-18859

**Singapore** ■ 65-67775366

**Spain** ■ 91-630-7050

**Sweden** ■ 020-790282

**Switzerland** ■ 055-254-22-11

**UK** ■ 01293-422-911

**USA** ■ 800-426-8157



---

Sample & Assay Technologies