HotStarTaq[®] DNA Polymerase and QIAGEN[®] Taq DNA Polymerase

For high PCR specificity with minimal PCR optimization

HotStarTaq[®] DNA Polymerase provides high specificity in hot-start PCR and QIAGEN[®] Taq DNA Polymerase provides optimal results in a wide range of PCR applications.

HotStarTaq DNA Polymerase offers:

- High PCR specificity due to the combination of enzyme with unique QIAGEN PCR Buffer
- Reduced nonspecific amplification provided by the stringent built-in hot start
- Minimal optimization saving time and money
- Easy handling simple reaction setup at room temperature

Increased PCR specificity

HotStarTaq DNA Polymerase, together with unique QIAGEN PCR Buffer, provides high specificity in hot-start PCR. Amplification of nonspecific products, primer–dimer formation, and background are minimized in every PCR cycle (Figure 1).

HotStarTaq DNA Polymerase

HotStarTaq DNA Polymerase, a modified form of QIAGEN *Taq* DNA Polymerase, is supplied in an inactive state that has no polymerase activity at ambient temperatures. This prevents extension of nonspecifically annealed primers and primer-dimers formed at low temperatures during PCR setup and the initial PCR cycle. HotStarTaq DNA Polymerase is activated by a 15-minute incubation at 95°C that can be incorporated into any existing thermal-cycler program. The stringent hot start provided by HotStarTaq DNA Polymerase in combination with unique QIAGEN PCR Buffer results in the highest PCR specificity compared with other hot-start enzymes (Figure 2).









Figure 2. Different hot-start methods. A 497 bp fragment was amplified from 50 copies of an HIV-*pol*-gene construct which had been added to 1 µg human genomic DNA. Different hot-start enzymes were employed: HotStarTaq DNA Polymerase from QIAGEN; hot-start enzymes from Suppliers A_{II} and R; *Taq*-antibody mixture from Supplier I; or an enzyme with no hot start. Arrow indicates the specific PCR product. Equal volumes of the reaction were analyzed on a 2% agarose gel. **M**: markers.

Minimal PCR optimization with unique PCR Buffer

QIAGEN PCR Buffer, supplied with both HotStarTaq DNA Polymerase and *Taq* DNA Polymerase, has been developed to eliminate the need for optimization of individual primer-template systems, saving time and money. The balanced combination of KCl and $(NH_4)_2SO_4$ in the buffer promotes specific primer-template annealing. Simultaneously, nonspecific annealing is reduced, maximizing yields of specific PCR product. Specific amplification is maintained over a wide range of temperatures and Mg²⁺ concentrations, without the need for timeconsuming optimization (Figure 3).

Optimal results with QIAGEN Taq DNA Polymerase

Recombinant QIAGEN *Taq* DNA Polymerase in combination with the unique PCR Buffer has been successfully used in a variety of applications, including standard amplification, differential display, and RT-PCR.

Q-Solution® for amplification of difficult templates

HotStarTaq DNA Polymerase and *Taq* DNA Polymerase are provided with Q-Solution — an innovative PCR additive. Q-Solution facilitates the amplification of templates that are GC-rich or have extensive secondary structures by modifying the melting behavior of DNA. Use of this unique reagent will often enable or improve PCR of difficult sequences (Figure 5).

Table 1. Features and benefits of QIAGEN enzymes for PCR*

Enzyme	Hot start for high specificity	Unique buffer system for minimal optimization	Amplification of GC-rich templates
HotStarTaq DNA Polymerase	1	\checkmark	\checkmark
Taq DNA Polymerase	-	1	1

* ProofStart® DNA Polymerase provides high-fidelity amplification. Visit **www.qiagen.com** for further information on high-fidelity PCR.

Figure 3. Tolerance to wide annealing temperatures and variable Mg²⁺ concentrations. PCR amplification at the indicated annealing temperatures **A** and Mg²⁺ concentrations **B** using QIAGEN PCR Buffer and QIAGEN *Taq* DNA Polymerase (similar results are obtained using HotStarTaq DNA Polymerase). The same PCR was performed in parallel using PCR buffer and *Taq* DNA polymerase from another supplier (**Supplier A₁₁**). **A** Amplification of the single-copy human cystic fibrosis gene and **B** the single-copy human prion protein gene. **M**: markers.

Convenient master-mix format

The HotStarTaq Master Mix Kit and Taq PCR Master Mix Kit provide:

- Easy reaction setup convenient master-mix format
- Less pipetting with minimal risk of contamination
- Time-savings eliminates tedious handling steps
- Minimal optimization due to unique QIAGEN PCR buffer

Fast, easy, and convenient PCR setup

With the HotStarTaq Master Mix Kit and the *Taq* PCR Master Mix Kit, there is no need for separate pipetting of individual components, minimizing the risk of contamination. Lengthy calculations are avoided and setting up amplification reactions is fast and easy (Figure 6). Since errors due to pipetting variabilities and miscalculation are reduced, PCR results are highly reproducible. In addition, the hot start provided by HotStarTaq DNA Polymerase in the HotStarTaq Master Mix Kit leads to increased PCR specificity and amplification reactions can be set up at room temperature. To choose which enzyme is right for your application, see Table 1. Specifications for HotStarTaq DNA Polymerase and *Taq* DNA Polymerase are given in Table 2.







Table 2. Specifications for *Taq* DNA Polymerase and HotStarTaq DNA Polymerase

Concentration:	5 units/µl	
Recombinant enzyme:	Yes	
Substrate analogues:	dNTP, ddNTP, dUTP, biotin-11-dUTP, DIG-11-dUTP, fluorescent dNTP/ddNTP	
Extension rate:	2–4 kb/min at 72°C	
Half-life:	10 min at 97°C 60 min at 94°C	
Amplification efficiency:	≥10 ⁵ fold	
5'→3' exonuclease activity:	Yes	
Extra A addition:	Yes	
3'→5' exonuclease activity:	No	
Contaminating nucleases:	No	
Contaminating RNases:	No	
Self-priming activity:	No	

Figure 6. Comparison of standard PCR setup, and PCR setup using *Taq* PCR Master Mix and HotStarTaq Master Mix from QIAGEN.

Ordering Information

Product	Contents	Cat. no.
HotStarTaq DNA Polymerase – for high PCR	specificity	
HotStarTaq DNA Polymerase* (250 U)	250 units HotStarTaq DNA Polymerase, 10x PCR Buffer, [†] 5x Q-Solution, 25 mM MgCl ₂	203203
HotStarTaq DNA Polymerase* (1000 U)	4 x 250 units HotStarTaq DNA Polymerase, 10x PCR Buffer, [†] 5x Q-Solution, 25 mM MgCl ₂	203205
HotStarTaq Master Mix Kit – premixed solut	ion for high PCR specificity	
HotStarTaq Master Mix Kit (250 U)	3 x 0.85 ml HotStarTaq Master Mix [‡] containing 250 units HotStarTaq DNA Polymerase total, 2 x 1.7 ml distilled water	203443
Taq DNA Polymerase – for standard and sp	ecialized PCR applications	
Taq DNA Polymerase (250 U)*	250 units Taq DNA Polymerase, 10x PCR Buffer,† 5x Q-Solution, 25 mM MgCl ₂	201203
Taq DNA Polymerase (1000 U)	1000 units <i>Taq</i> DNA Polymerase, 10x PCR Buffer,† 5x Q-Solution, 25 mM MgCl ₂	201205
Taq PCR Core Kit (250 U)	250 units <i>Taq</i> DNA Polymerase, 10x PCR Buffer,† 5x Q-Solution, 25 mM MgCl ₂ , dNTP Mix§	201223
Taq PCR Master Mix Kit — premixed solutio	n for covenient PCR setup	
Taq PCR Master Mix Kit (250 U)	3 x 1.7 ml <i>Taq</i> PCR Master Mix [‡] containing 250 units <i>Taq</i> DNA Polymerase total, 3 x 1.7 ml distilled water	201443

* Kits containing up to 25,000 units HotStarTaq DNA Polymerase or Taq DNA Polymerase are available; please inquire.

[†] Contains 15 mM MgCl₂

 ‡ Provides a final concentration of 1.5 mM MgCl_ and 200 μM each dNTP

 $\ensuremath{\$}$ Contains 10 mM each dNTP

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at **www.qiagen.com** or can be requested from QIAGEN Technical Services or your local distributor.

Visit www.qiagen.com/goto/PCR or contact QIAGEN today to discover how easy and specific your PCR can be!

Trademarks: QIAGEN®, Sample to Insight[®], HotStarTaq[®], ProofStart[®], Qa-Solution[®] (QIAGEN Group). Purchase of QIAGEN products for PCR is accompanied by a limited license to use them in the Polymerase Chain Reaction (PCR) process for research and development activities in conjunction with a thermal cycler whose use in the automated performance of the PCR process is covered by the up-front license fee, either by payment to Applied Biosystems or as purchased, i.e. an authorized thermal cycler. The PCR process is covered by U.S. Patents 4,683,195 and 4,683,202 and foreign equivalents owned by Hoffmann-La Roche AG. © 2015 QIAGEN, all rights reserved. PROM-9014-001

Ordering www.qiagen.com/contact | Technical Support support.qiagen.com | Website www.qiagen.com