

Guidelines for installation of Q Protocols for QIAcube[®] HT Kits on the QIAcube HT and the QIAxtractor[®]


These guidelines outline the procedure for the installation of QIAcube HT Q Protocols onto your computer. When the QIAcube HT Software is installed, double-click on the Q Protocol MSI-Windows[®] Installer Package to identify the installation path. The Setup Wizard uses this path to install the Q Protocols in the correct folder. QIAcube HT Q Protocols are automatically saved in the "QProtocols" folder of the QIAcube HT Software (C:\Program Files\QIAcubeHT\QProtocols).

Installing Q Protocols

1. Go to www.qiagen.com/p/QIAcubeHT
2. Select the "Resources" tab.
3. Under the "Protocol Files" header, double-click on the Q Protocol to be installed.
Note: For more information about a Q Protocol, open the corresponding "Q Protocol Note" in the folder/header "Protocol Files".
4. Save the Q Protocol MSI Installation Package on any drive of your computer.
5. Double-click on the Q Protocol MSI Installation Package (e.g., QIAamp[®] 96 DNA QIAcube HT V1.msi).
6. If the QIAcube HT Software is installed on the computer, installation of the selected Q Protocol will start.
Note: For information on how to install QIAcube HT software, please refer to the *QIAcube HT User Manual* which can be downloaded on www.qiagen.com/p/QIAcubeHT, under the "Resources" tab.
7. Click "Finish" to complete and exit the Setup Wizard.



8. The Q Protocol has now been installed in the QProtocols folder of the QIAcube HT Software.

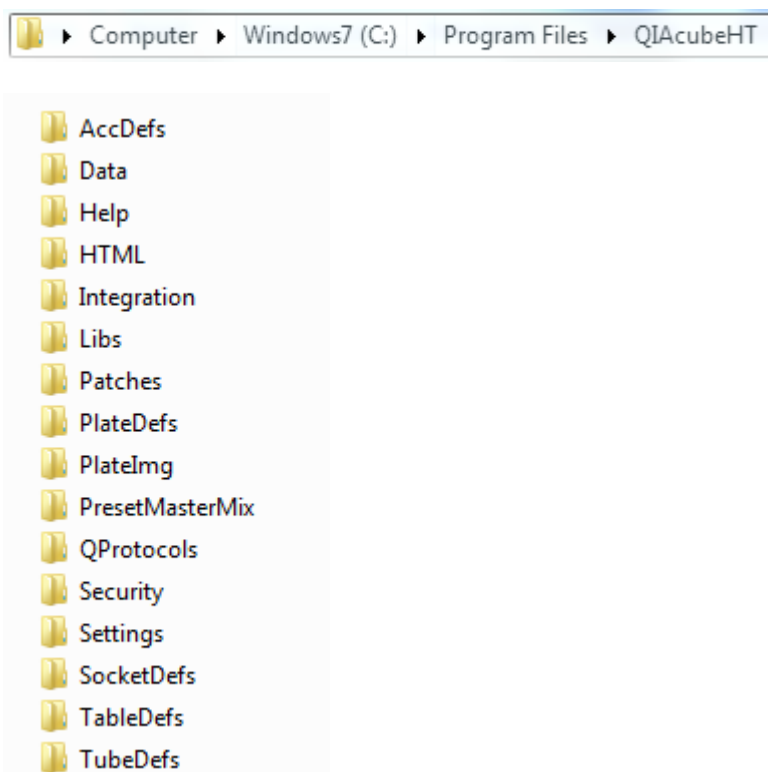

Note: For details on how to use the Q protocols, please refer to the *QIAcube HT User Manual*, which can be downloaded on www.qiagen.com/p/QIAcubeHT, under the "Resources" tab.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or use manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, QIAamp®, QIAcube®, QIAxtractor® (QIAGEN Group); Windows® (Microsoft Corporation).

June-13 © 2013 QIAGEN, all rights reserved.

www.qiagen.com

Australia ■ 1-800-243-800

Austria ■ 0800/281010

Belgium ■ 0800-79612

Canada ■ 800-572-9613

China ■ 021-51345678

Denmark ■ 80-885945

Finland ■ 0800-914416

France ■ 01-60-920-930

Germany ■ 02103-29-12000

Hong Kong ■ 800 933 965

Ireland ■ 1800 555 049

Italy ■ 800 787980

Japan ■ 03-5547-0811

Korea (South) ■ 1544 7145

Luxembourg ■ 8002 2076

The Netherlands ■ 0800 0229592

Norway ■ 800-18859

Singapore ■ 65-67775366

Spain ■ 91-630-7050

Sweden ■ 020-790282

Switzerland ■ 055-254-22-11

UK ■ 01293-422-911

USA ■ 800-426-8157



Sample & Assay Technologies