



July 2025

Important Note

EZ1&2® ccfDNA Kit (48) (cat. no. 954854)

Dear valued EZ1&2 ccfDNA Kit customer,

We would like to inform you that we have identified a bug in EZ2 Standard Protocol Package version 10 and 11 in the following experiment setup:

Parameter	Selection value
Application	DNA
Kit	ccfDNA Kit
Protocol	ccfDNA from plasma or serum
Sample volume	10 mL
Elution volume	75 µL

This bug results in an elution volume lowered by approximately 25 µL, but it will not have an impact on the quality or yield. Other settings and protocols are not affected.

At QIAGEN, we strive to provide the highest level of performance with our instruments and protocols. Therefore, the fix for this bug will be available with the upcoming Protocol Package version 12 which can be downloaded from the EZ2 Connect website in August 2025. For early access to the updated protocol, please contact your local QIAGEN representative or our Technical Support Center at support.qiagen.com

Sincerely,

Your EZ2 Connect Team

Trademarks: QIAGEN®, Sample to Insight®, EZ1&2® (QIAGEN Group). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law. 07/2025 HB-3730-001 © 2025 QIAGEN, all rights reserved.