

PyroMark® Q24 Vacuum Workstation quick-start guide


This Technical Information summarizes the immobilization and preparation of PCR products for Pyrosequencing® on the PyroMark Q24 Instrument. Before using the PyroMark Q24 Vacuum Workstation, carefully read Section 5.3 of the *PyroMark Q24 User Manual* and pay particular attention to the safety information.

Immobilizing the PCR products

1. Make a master mix according to the flowchart to the right.

Note: Before pipetting, gently shake the bottle of streptavidin-coated Sepharose® beads* to ensure a homogenous suspension.


2. Depending on the sample volume, dispense 60–75 µl master mix into each necessary well of a PCR plate to give a total volume of 80 µl per well.
3. Add 5–20 µl PCR product to each well.
4. Seal the wells with strip caps and agitate the PCR plate at 1400 rpm for 5–10 min at room temperature (15–25°C) using an orbital shaker.

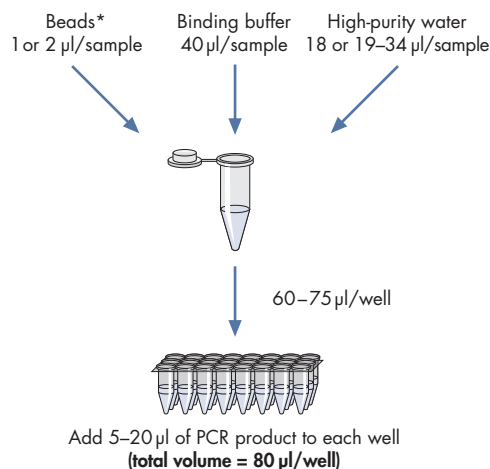

Preparing the master mix to immobilize the PCR products.

Separation of DNA strands and release of samples into the PyroMark Q24 Plate

1. Dilute the sequencing primers to 0.3 µM with PyroMark Annealing Buffer (cat. no. 979009), and dispense 25 µl into each necessary well of a PyroMark Q24 Plate. Position the plate on the workstation.
2. Fill the workstation troughs according to the diagram to the right.
3. Start the pump and apply vacuum to the tool by opening the switch. Flush the filter probes with high-purity water (Milli-Q® 18.2 MΩ x cm or equivalent) in trough 5. Refill the trough with fresh high-purity water for use in step 12.
4. Position the PCR plate on the workstation. Ensure that both plates are in the same orientation as when the samples were loaded.
5. With the vacuum switch ON, lower the vacuum tool into the wells of the PCR plate for 15 s to capture the beads with PCR product.

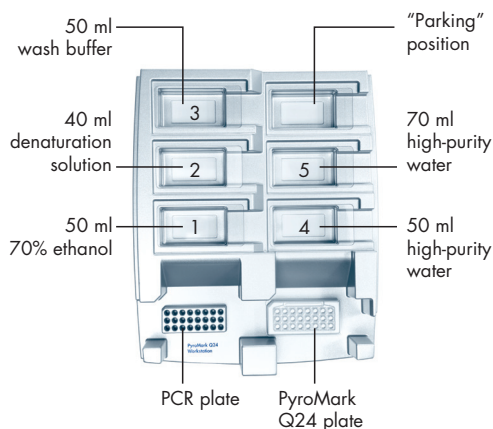

Filling the workstation troughs.

* Streptavidin Sepharose High Performance (34 µm, 5 ml, GE Healthcare). Check the lot number of the Streptavidin Sepharose High Performance. For lot number 10057037 or higher, use 1 µl in the master mix. For lot numbers lower than 10057037, use 2 µl. This is not a complete list of suppliers and does not include many important vendors of biological supplies.





CAUTION:

PyroMark Denaturation Solution contains sodium hydroxide, which is an eye and skin irritant.

Always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, see the *PyroMark Q24 User Manual* (and the MSDS).



Vacuum switch ON.



Vacuum switch OFF.

6. With the vacuum ON, flush the tool with 70% ethanol (trough 1) for 5 s.
7. With vacuum ON, flush the tool with denaturation solution (trough 2) for 5 s.
8. With vacuum ON, flush the tool with wash buffer (trough 3) for 10 s.
9. With vacuum ON, raise the vacuum tool to beyond 90° vertical for 5 s.
10. Align the vacuum tool with the PyroMark Q24 Plate and switch the vacuum OFF. Lower the vacuum tool into the wells and gently shake from side to side to release the beads.
11. With the vacuum OFF, agitate the vacuum tool in high-purity water (trough 4) for 10 s.
12. Switch the vacuum ON and flush the filter probes with high-purity water (trough 5) for 5 s.
13. Raise the vacuum tool to beyond 90° vertical for 5 s, then switch the vacuum OFF and store the tool in the "Parking" position.

Annealing sequencing primers to DNA strands

Place the PyroMark Q24 Plate in a prewarmed PyroMark Q24 Plate Holder. Heat the Pyrosequencing samples on a heating block at 80°C for 2 minutes.

Remove the plate from the holder and allow the samples to cool to room temperature (15–25°C) for at least 5 minutes. The cooled plate can now be processed in the PyroMark Q24.

Cleaning the vacuum workstation

Liquid waste and solutions remaining in the troughs of the vacuum workstation should be appropriately discarded at the end of the day. For details, see the *PyroMark Q24 User Manual*.

For up-to-date licensing information and product-specific disclaimers, see the respective *PyroMark Q24 User Manual*. The *PyroMark Q24 User Manual* is available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN®, PyroMark®, Pyrosequencing® (QIAGEN Group); Milli-Q® (Millipore Corporation); Sepharose® (GE Healthcare).
1074528 12/2012 © 2012 QIAGEN, all rights reserved.

Australia = 1-800-243-800
Austria = 0800-281011
Belgium = 0800-79612
Brazil = 0800-557779
Canada = 800-572-9613
China = 800-988-0325
Denmark = 80-885945

Finland = 0800-914416
France = 01-60-920-930
Germany = 02103-29-12000
Hong Kong = 800 933 965
India = 1-800-102-4114
Ireland = 1800 555 049
Italy = 800-787980

Japan = 03-6890-7300
Korea (South) = 080-000-7145
Luxembourg = 8002 2076
Mexico = 01-800-7742-436
The Netherlands = 0800-0229592
Norway = 800-18859
Singapore = 1800-742-4368

Spain = 91-630-7050
Sweden = 020-790282
Switzerland = 055-254-22-11
Taiwan = 0080-665-1947
UK = 0808-234-3665
USA = 800-426-8157
www.qiagen.com

