

# EZ1<sup>®</sup> Virus Mini Kit v2.0

The EZ1 Virus Mini Kit v2.0 (cat. no. 955134) can be stored at room temperature (15–25°C) until the indicated expiration date. Do not freeze the reagent cartridges.

## Further information

- *EZ1 Virus Handbook*: [www.qiagen.com/HB-0113](http://www.qiagen.com/HB-0113)
- Safety Data Sheets: [www.qiagen.com/safety](http://www.qiagen.com/safety)
- Technical assistance: [support.qiagen.com](http://support.qiagen.com)

## Notes before starting

- The EZ1 Virus Mini Kit v2.0 requires use of the EZ1 Advanced XL Virus Card v2.0 with the EZ1 Advanced XL, or use of the EZ1 Advanced Virus Card v2.0 with the EZ1 Advanced, or use of the EZ1 Virus Card v2.0 with the BioRobot<sup>®</sup> EZ1.
- The buffer in well 1 of the reagent cartridge (i.e., the well that is nearest to the front of the EZ1 instrument when the reagent cartridge is loaded) may form a precipitate upon storage. If necessary, redissolve by warming at 37°C and then place at room temperature (15–25°C).
- EZ1 instruments should only be switched on after an EZ1 Card is inserted. Make sure that the EZ1 Card is completely inserted, otherwise essential instrument data could be lost, leading to a memory error. EZ1 Cards should not be exchanged while the instrument is switched on.
- Before loading reagent cartridges into the EZ1 instrument, invert the cartridges three times to mix the magnetic particles then tap to deposit the reagents at the bottom of their wells. Check that the magnetic particles are completely resuspended.

1. If necessary, pretreat samples according to the appropriate procedure, as described in the *EZ1 Virus Handbook*. See Table 1 for information on which sample types require pretreatment as well as accessories required for pretreatment.
2. Carry out nucleic acid purification using an EZ1 instrument with the appropriate EZ1 Card, according to the *EZ1 Virus Handbook*.

**Note:** Specific guidance regarding the amount of starting material for different sample types is provided in the *EZ1 Virus Handbook*.

**Table 1. Sample materials covered by the EZ1 Virus Mini Kit v2.0**

Sample type	Pretreatment needed	Pretreatment accessories	Ordering information
Serum	No	–	–
Plasma	No	–	–
Cerebrospinal fluid	No	–	–
Urine*	Yes*	Buffer ATL	Cat. no. 939011
Whole Blood*	Yes*	Buffer ATL	Cat. no. 939011
Stool*	Yes*	Buffer ASL	Cat. no. 19082
Transport media	No	–	–
Dried swabs*	Yes*	Buffer ATL	Cat. no. 939011
Respiratory samples*	Yes*	Sputasol	<b>www.oxid.com</b>

\* Information about specific pretreatment procedures is provided in the *EZ1 Virus Handbook*.



Scan QR code for handbook.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.

Trademarks: QIAGEN®, Sample to Insight®, BioRobot®, EZ1® (QIAGEN Group). 1101295 03/2016 HB-0787-002 © 2016 QIAGEN, all rights reserved.