

March 2011

RT² miRNA qPCR Assay Handbook

For miRNA expression analysis



Sample & Assay Technologies

QIAGEN Sample and Assay Technologies

QIAGEN is the leading provider of innovative sample and assay technologies, enabling the isolation and detection of contents of any biological sample. Our advanced, high-quality products and services ensure success from sample to result.

QIAGEN sets standards in:

- Purification of DNA, RNA, and proteins
- Nucleic acid and protein assays
- microRNA research and RNAi
- Automation of sample and assay technologies

Our mission is to enable you to achieve outstanding success and breakthroughs. For more information, visit www.qiagen.com.

Product Use Limitations

RT² miRNA qPCR Assays are intended for molecular biology applications. This product is not intended for the diagnosis, prevention, or treatment of a disease.

All due care and attention should be exercised in the handling of the products. We recommend all users of QIAGEN products to adhere to the NIH guidelines that have been developed for recombinant DNA experiments, or to other applicable guidelines.

CONTENTS

I.	Background and Introduction	4
II.	Materials Provided	6
III.	Additional Materials Required	6
IV.	Protocol	7
	A. RNA Preparation and Quality Control	9
	B. RT ² miRNA First Strand Kit	11
	C. Performing Real-Time PCR	12
	D. Data Analysis: $\Delta\Delta C_t$ Method	15
	E. Data Analysis: Optional Standard Curve Method	17
VI.	Troubleshooting and Frequently Asked Questions	18

I. Background and Introduction

MicroRNA (miRNA), first discovered in *C. elegans* in early 1990s, are a class of naturally occurring small RNA molecules, generally 17 to 30 nucleotides long, processed from much larger stem-loop structures (pri-miRNA and pre-miRNA). Mature miRNA is recognized by and guides the RNA-Induced Silencing Complex (RISC) to specific messenger RNA transcript sequences, causing either translation repression or mRNA degradation. The expression and function of miRNA brings another layer of gene expression regulation to an already complex network. Each miRNA can regulate multiple mRNA targets, and a single target mRNA may be regulated by multiple miRNA sequences. Researchers are currently correlating miRNA expression profiles to biological phenotypes in attempts to better understand miRNA-based gene regulation. Real-time, reverse transcription PCR, or qRT-PCR, is the most sensitive and reliable method for nucleic acid expression analysis currently available. However, the small size and high degree of similarity among miRNA sequences makes quantitative analyses of their abundance very challenging.

The RT² miRNA qPCR Assays take advantage of primer design and assay formulation technologies to analyze the expression of any mature miRNA sequence in the human, mouse, or rat microRNA genome (miRNome), as annotated by the Sanger miRBase Release 14, by SYBR[®] Green based real-time RT-PCR. The assays rely on combining poly(A) tailing and a universal reverse transcription in one first strand cDNA synthesis reaction. Our proprietary primer design algorithm also distinguishes miRNA family members with single nucleotide mismatches. Each assay is experimentally validated with an *in vitro* assay to insure the amplification of a single product with at least consistently high efficiency under standardized cycling conditions. The high performance of every RT² miRNA qPCR Assay is certified and 100% guaranteed when performed using the complete miRNA PCR Assay System that includes the RT² miRNA First Strand Kit and RT² SYBR Green qPCR Mastermix.

To complete the miRNA PCR Assay procedure, reverse transcribe your experimental small RNA samples into first strand cDNA, the template for the PCR, using our RT² miRNA First Strand Kit. (See Figure 1 for an overview of the miRNA PCR Assay procedure.) Then, mix the template with one of our instrument-specific and ready-to-use RT² SYBR Green qPCR Mastermixes. Set up your assays, perform PCR, and finally determine relative expression with your real-time instrument and the $\Delta\Delta C_t$ method. The simplicity of the RT² miRNA qPCR Assays makes them accessible for routine use in every research laboratory.


Benefits of the RT² miRNA qPCR Assays:

- **High Performance:** Each assay is experimentally validated to amplify a single amplicon with uniform PCR efficiency.
- **Complete Genome Coverage:** Assays are available for over 400 majority miRNA sequences. The uniform PCR efficiency and conditions allow easy, scalable transitions from analyzing single to multiple miRNA sequences.
- **Simple & Convenient:** Universal reverse transcription converts all miRNA into cDNA in a one-step reaction. Assays deliver guaranteed performance when used with


optimized instrument-specific RT² SYBR Green qPCR Mastermixes for ABI, Bio-Rad[®], Stratagene[®] and other real-time PCR instruments.

How miRNA PCR ARRAYS WORK

1. Convert miRNA to cDNA.

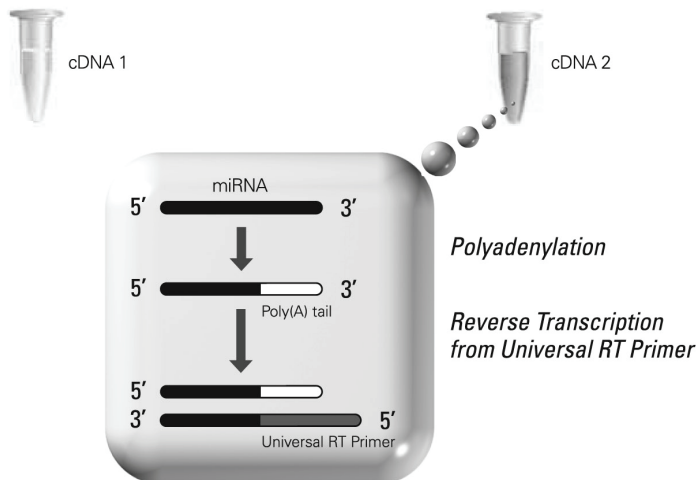

2. Add cDNA to RT² qPCR Master Mix. Aliquot Mixture Across PCR Array.

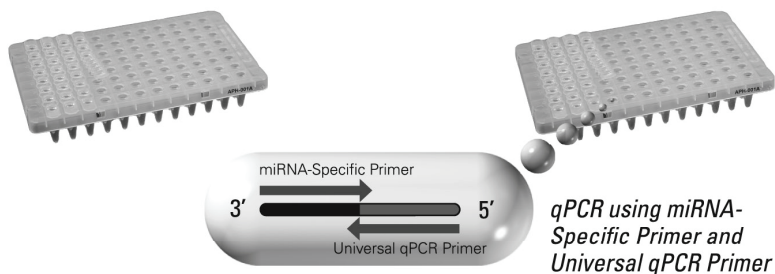

3. Run in Your Real-Time PCR Instrument.



4. Data Analysis.

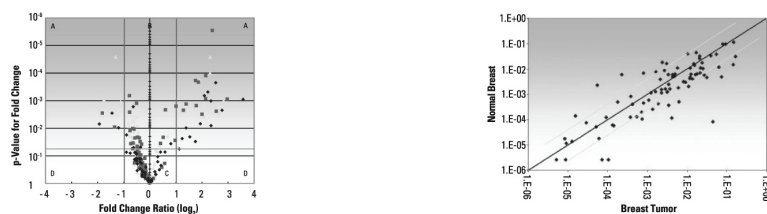


Figure 1: Overview of the RT² miRNA qPCR Assay procedure.

II. Materials Provided:

RT² miRNA qPCR Assay, one tube, 200-reaction scale

- Primer pair includes one miRNA-specific primer and one universal primer.
- One Product Specification Sheet

Storage Conditions:

Primer Sets are shipped at ambient temperature but must be stored at –20°C where they are guaranteed as is for 6 months from the date received.

III. Additional Materials Required:

A. miRNeasy Mini Kit (Cat. No. 217004)

B. RT² miRNA First Strand Kit (12) (Cat. No. 331401)

MANDATORY for a Complete and Successful Experiment

The universal primer in each assay is only compatible with the sequence added to the cDNA template by the primer in this kit.

C. RT² SYBR Green qPCR Mastermix

MANDATORY for a Complete and Successful Experiment

Be sure to pick the correct one for the instrumentation in your laboratory.

RT² SYBR Green ROX™ qPCR Mastermix:

Specifically designed for: All ABI and Stratagene Instrumentation and Eppendorf® Mastercycler® ep *realplex* Instruments with ROX filter set

Catalog Number Size

330520 For 200 25-μL reactions (2.5 mL)

330529 For 2000 25-μL reactions (25.0 mL)

RT² SYBR Green Fluor qPCR Mastermix:

Specifically designed for BioRad iCycler®, MyiQ™, and iQ™5 Instrumentation

Catalog Number Size

330510 For 200 25-μL reactions (2.5 mL)

330519 For 2000 25-μL reactions (25.0 mL)

RT² SYBR Green qPCR Mastermix:

Specifically designed for instrumentation that does not require a reference dye:

BioRad (MJ Research) Opticon, Opticon 2, and Chromo™ 4

Roche® LightCycler® 480 System

Eppendorf Mastercycler ep *realplex* Instruments without ROX filter set

Catalog Number Size

330500 For 200 25-μL reactions (2.5 mL)

330509 For 2000 25-μL reactions (25.0 mL)

D. Equipment:

1. For recommendations on specific real-time instrumentation (thermal cyclers with fluorescent detection), see the list of master mixes and plate formats above.
2. Calibrated Pipettors
3. RNase / DNase-free pipette tips and tubes

IV. Protocol:

Please read through this entire protocol before beginning your experiment. RNA samples are very sensitive to RNase digestion; therefore, wear gloves and maintain an RNase-free work area while performing this protocol.

NOTE: Master Mix and Reverse Transcription Kit Considerations

The performance of our RT² miRNA qPCR Assays is only guaranteed with our RT² SYBR Green qPCR Mastermixes and the RT² miRNA First Strand Kit. Therefore, the use of the complete miRNA qPCR Assay System is absolutely essential for obtaining any real-time PCR profiling results.

The chemically-modified and tightly controlled HotStart enzyme and other proprietary chemical components in our RT² SYBR Green qPCR Mastermixes provide more accurate SYBR Green results by preventing the amplification of primer dimers and other non-specific products. They also help insure high amplification efficiencies. When we test other sources of enzymes with our RT² miRNA qPCR Assays, we frequently see primer dimers and other non-specific products that confound SYBR-Green based real-time PCR detection. Because each instrument uses a different reference dye to normalize their optics, be sure that you use the correct master mix for the instrumentation in your laboratory.

The universal primer in every assay of the RT² miRNA qPCR Assays is specific only for the unique and proprietary sequence incorporated into the cDNA by the universal RT primer in the RT² miRNA First Strand Kit. The RT² miRNA qPCR Assays cannot detect cDNA generated from miRNA using other sources of miRNA first strand synthesis or miRNA reverse transcription kits. Similarly, cDNA generated from the RT² miRNA First Strand Kit cannot be used with other sources of real-time qPCR assays for miRNA.

The RT² miRNA First Strand Kit also includes a proprietary buffer to preferentially reverse transcribe miRNA into cDNA. The buffer components and the magnesium concentration in our RT² miRNA First Strand Kit are also more compatible with our PCR master mixes than other enzymes or kits providing the RT² miRNA qPCR Assays with maximum levels of sensitivity with ng amounts of small RNA.

NOTE: Preparing a Workspace Free of DNA Contamination

For accurate and reproducible PCR assay results, it is very important to avoid contamination of the assay with foreign DNA. Any DNA contamination will artificially inflate the SYBR Green signal yielding skewed expression profiles and false positive signals. The most common sources of DNA contamination are the products of previous experiments spread into the air of your working environment. Please follow the recommendations below on how to set up and maintain a working environment free of DNA contamination.

1. Wear gloves throughout the procedure. Use only fresh PCR-grade reagents (H₂O) and lab ware (tips and tubes).
2. Physically separate the workspaces used for PCR setup and post-PCR processing or non-PCR operations. Decontaminate your PCR workspace and lab ware (pipettor barrels, tube racks, etc.) before each new use with UV light to render any contaminating DNA ineffective in PCR through the formation of thymidine dimers or with 10% bleach to chemically inactivate and degrade any DNA.
3. Close all tubes containing PCR products once you are finished adding or removing volumes. Before discarding any lab ware (tips or tubes) containing PCR products or other DNA, treat with 10% bleach.
4. Do not open any previously run and stored PCR plate. Removing the thin-wall 8-cap strips or the adhesive film from PCR plates releases PCR product DNA into the air where it will contaminate and confound the results of future real-time PCR experiments.

A. RNA Preparation and Quality Control:

High quality RNA is **ESSENTIAL** for obtaining good real-time PCR results.

The most important prerequisite for any miRNA expression analysis experiment is consistent, high-quality small RNA from every experimental sample. Therefore, the sample handling and RNA isolation procedures are critical to the success of the experiment. Residual traces of proteins, salts or other contaminants will either degrade the RNA or decrease the efficiency of (if not block completely) the enzyme activities necessary for optimal reverse transcription and real-time PCR performance.

Although other real-time PCR-based miRNA detection methods call for the use of total RNA that contains miRNA as the starting input material, we highly recommend starting with small RNA (< 200 nucleotides) to reduce background noise for optimal sensitivity.

1. Recommended RNA Preparation Methods:

High quality small RNA for your real-time PCR experiment must be prepared using one of the following methods, each specific for your biological sample:

Cultured Cells OR Tissue Samples: Use the miRNeasy Mini Kit (cat. no. 217004).

- a. Extract RNA from the tissue using the the QIAzol[®] Lysis Reagent included in the miRNeasy Mini Kit. Be sure to use a sufficient amount of QIAzol Lysis Reagent. During homogenization, add a volume of reagent at least ten times greater than the tissue volume.
- b. Then further clean-up and enrich the small RNA following the protocol provided in the miRNeasy Mini Handbook.

Whole Blood Samples:

Remove red blood cells (RBC) from whole blood samples using a density gradient centrifugation medium (for example, Lymphoprep, Greiner Bio-One, Catalog # 1031966). Isolate small RNA from the white blood cell fraction as described for cultured cells above.

Previously Isolated Total RNA:

If you have already prepared total RNA from any biological source material, be sure that it used a phenol-based method (such as TRIzol, RNeasy, etc.). Then, enrich for small RNA using the miRNeasy Mini Kit (cat. no. 217004).

For Other Biological Samples: Refer to the existing literature to find isolation protocols for high-quality RNA from other biological samples or contact one of our Technical Support representatives.

For best results from the PCR Assays, all RNA samples should be suspended in the RNase-free water provided with the RNA Isolation kit, or alternatively in RNase-free 10 mM Tris buffer pH 8.0. DO NOT use DEPC-treated water!

2. RNA Quality Control:

For best results from the PCR Assays, all RNA samples should also demonstrate consistent quality according to the following criteria:

a. RNA Concentration and Purity by UV Spectrophotometry

NOTE: Prepare dilutions and measure absorbance in 10 mM Tris, pH 8.0 buffer. The spectral properties of nucleic acids are highly dependent on pH.

- i) $A_{260}:A_{230}$ ratio should be greater than 1.7.
- ii) $A_{260}:A_{280}$ ratio should be greater than 2.0.
- iii) Concentration by A_{260} should be greater than 1 μg / ml small RNA

b. RNA Quality & Integrity

Characterize ~10 ng of the small RNA on an Agilent® Bioanalyzer using an RNA 6000 Nano LabChip®. The RNA should contain a single sharp peak at a low molecular weight with no smearing and no additional peaks at higher molecular weights.

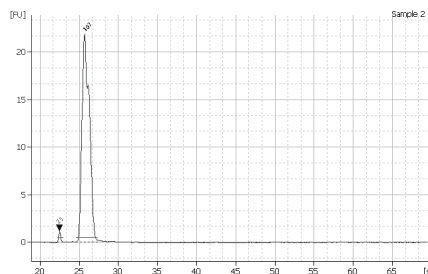


Figure 3: Good Small RNA Band Integrity Is Important for Best Results from the miRNA qPCR Array. Displayed is an Agilent BioAnalyzer electropherogram of a high-quality small RNA preparation showing a sharp low molecular weight peak without a shoulder.

3. Amount Considerations:

The RT² miRNA qPCR Assay System yields relative miRNA expression profiles with as little as 1 ng or as much as 200 ng of small RNA. However, the optimal amount of starting material depends upon the relative abundance of the sequences of interest. Lower abundance sequences require more RNA; higher abundance sequences require less RNA. Greater amounts of input RNA yield a greater number of positive calls; that is, sequences expressed in the linear dynamic range of the method. Lower amounts of input RNA yield a smaller number of positive calls and increase false negative calls.

The use of the RT² miRNA First Strand Kit maximizes the number of positive calls at low amounts of RNA over other sources of miRNA reverse transcriptase and first strand synthesis kits. For successful results and maximum positive call rates, we recommend that first time users try starting with 100 ng of small RNA. It is also important to use a consistent amount of RNA for each sample to be characterized and compared.

B. RT² miRNA First Strand Kit

NOTE: RNA samples must meet the standards of integrity and purity from protein and organics contamination described on the previous two pages.

NOTE: The buffer components of the reverse transcription reaction using this kit are more compatible with our PCR master mixes than other sources of kits for first strand cDNA synthesis kits or reverse transcriptases.

NOTE: Thaw 5x miRNA RT Buffer 2 on ice, briefly vortex the tube to mix the contents well, and centrifuge to the bottom of the tube before each use.

For each RNA sample, combine the following in a sterile PCR tube:

Small RNA	50 ng to 200 ng (or 0.5 – 2 µg total RNA)
miRNA RT Primer & ERC Mix	1.0 µL
5x miRNA RT Buffer 2	2.0 µL
miRNA RT Enzyme Mix	1.0 µL
miRNA Nucleotide Mix	1.0 µL
<hr/>	
RNase-free H ₂ O to a final volume of	10.0 µL

NOTE: Use the same amount of RNA in this reaction for every sample. First time users are recommended to start with 100 ng of small RNA. Do not use more than 400 ng of small RNA per reaction.

- b. Mix the contents gently with a pipettor followed by brief centrifugation.
- c. Incubate at 37°C for 2 hours.
- d. Heat at 95°C for 5 minutes to degrade the RNA and to inactivate the reverse transcriptase.
- e. Chill on ice for at least one minute and add 90 µl of RNase-free H₂O to each 10-µl of cDNA synthesis reaction. Mix well.
- f. Hold the finished First Strand cDNA Synthesis Reaction on ice until the next step or store overnight at -20°C.

C. Performing Real-Time PCR:

NOTE: *The use of RT² SYBR Green qPCR Mastermixes is critical for obtaining the most accurate results from the miRNA PCR Assays. Be sure to use the correct master mix for your instrument before continuing with this protocol (See Page 6).*

NOTE: *The accuracy and precision of your pipetting determines the consistency of your results. Be sure that all of your micro-pipettors are calibrated before beginning this procedure. Also, make sure to not introduce any bubbles into the PCR wells.*

NOTE: *If unsure of your RNA quality or isolation technique, examine the quality of your RNA before this step.*

1. Setup:

a. Experimental Samples

To insure the consistency of your results and that each experimental sample yields a reliably detectable C_t value, we recommend using undiluted template and a 1:10 dilution of your template in separate PCR assays. Also, prepare triplicate reactions for each template at each concentration.

For every experimental sample, prepare one set of reactions for every miRNA sequence of interest and for a single housekeeping sequence or a set of housekeeping sequences to normalize your raw data. Choose housekeeping sequence(s), such as small nuclear RNA species, known not to change their expression under your experimental conditions.

b. Positive and Negative Controls

Prepare a positive control reaction for each RT² miRNA qPCR Assay using cDNA template generated by reverse transcribing 2 µL of a 1 – 10 nM solution of synthetic RNA oligo with the same sequence(s) as the miRNA of interest.

To control for DNA contamination introduced during reaction setup, prepare a negative control reaction replacing template with water, the so-called No Template Control (NTC).

To monitor non-specific background amplification, perform one assay for each miRNA and each housekeeping sequence of interest using an equivalent volume of product from a reaction replacing the miRNA RT Enzyme Mix with RNase-free H₂O for each miRNA sample, the so-called No Reverse Transcription (NRT) control.

c. Optional Standard Curve Method

Generate one standard curve for each miRNA of interest. Also generate a standard curve for each housekeeping small nuclear RNA sequence to be used for data normalization.

To generate a standard curve, prepare a five-point series of five- or ten-fold dilutions in duplicate using cDNA template generated by reverse transcribing synthetic miRNA with the sequence(s) of interest.

2. Polymerase Chain Reactions

For each 25- μ l PCR, mix the following components in a PCR tube:

12.5 μ l	RT ² SYBR Green qPCR Mastermix, matched with your instrument*
10.5 μ l	ddH ₂ O
1.0 μ l	of either undiluted or diluted template
1.0 μ l	RT ² miRNA qPCR Assay primer set
<hr/>	
25.0 μ l	final volume

3. Performing Real-Time PCR Detection:

NOTE: *Be sure to follow the manufacturer's instructions for the proper operation and maintenance of your real-time instrument.*

- CAREFULLY** but tightly seal the PCR plate either with optical thin-wall 8-cap strips or optical adhesive film, according to the instrument manufacturer's instructions.

NOTE: *Be sure that no bubbles appear in any of the wells of the PCR plate. To remove bubbles, tap the plate gently on the bench top or centrifuge the plate briefly.*

- Place the plate on ice while setting up the PCR cycling program below.
- Place the plate in your real-time thermal cycler. If recommended by your instrument's user manual, use a compression pad on plates sealed with optical film.
- Enter and run the appropriate program for your real-time instrument (below). If prompted by your instrument software, select "Absolute Quantitation" to begin.

NOTE: *For additional help with instrument setup, see our Instrument-Specific Setup Instructions and Protocol Files at: www.sabiosciences.com/pcrarrayprotocolfiles.php*

Cycles	Duration	Temperature
1	10 minutes ¹	95°C
	15 seconds	95°C
40	30 to 40 seconds ^{2,3}	60°C
	30 seconds	72°C

¹ The 10-minute step at 95°C is required to activate the HotStart DNA polymerase.

² Detect and record SYBR Green fluorescence from every well during the annealing step of each cycle.

³ Different instruments need different lengths of time to detect the fluorescent signal. Choose the appropriate time for the annealing step (60°C) for your instrument.

- e. Calculate the threshold cycle (C_t) for each well using the instrument's software.
 - i. We highly recommend letting your instrument automatically define the Baseline values but manually setting the Threshold values.
 - ii. To define the Baseline, use the Linear View of the amplification plots and set the instrument to use the readings from cycle number two (2) through two (2) cycle values before the earliest visible amplification, usually around cycle number ten (10) but no more than 15.
 - iii. To define the Threshold Value, use the Log View of the amplification plots and place it above the background signal but within the lower half to one-third of the linear phase of the amplification plot.
 - iv. **IMPORTANT:** Ensure that the Thresholds are the same across all PCR runs in the same analysis.
 - v. Export the resulting threshold cycle values for all wells to a blank Excel spreadsheet for use with our Data Analysis Template Excel file.

4. Recommended Quality Control:

- a. Dissociation (Melting) Curve:

Run a melting curve program immediately after the above cycling program, and generate a first derivative dissociation curve for each well in the entire plate using your instrument's software. No more than one peak should appear in each reaction at temperatures greater than 70 °C. If your instrument does not have a default melting curve program, run the following program instead:

95°C, 1 min; 65°C, 2 min (OPTICS OFF); 65°C to 95°C at 2°C / min (OPTICS ON).

If you decide not to obtain the dissociation curve immediately, save the plates wrapped in aluminum foil at – 20°C as is, in case you need to perform this operation at a later point in time for troubleshooting purposes. When ready, simply warm the plate to room temperature, place it into your real-time instrument, and run the melting program described above.

NOTE: *Be sure to visually inspect the plate after the run for any signs of evaporation from any of the wells. If evaporation is observed, make a note of which wells so that you may qualify your data analysis appropriately.*

NOTE: *DO NOT open any previously run and stored PCR plate. Removing the thin-wall 8-cap strips or the adhesive film from PCR plates releases PCR product DNA into the air where it will contaminate and confound the results of future real-time PCR experiments. See also the Note on "Preparing a Workspace Free of DNA Contamination".*

D. Data Analysis: $\Delta\Delta C_t$ Method

1. Change all C_t values reported as greater than 35 or as N/A (not detected) to 35. At this point, any C_t value equal to 35 is considered a negative call.
2. Examine the threshold cycle values of the control wells.
 - a. Positive Control:
 - b. No Reverse Transcription (NRT):
A raw C_t value of 35 or greater indicates a lack of non-specific background.
A raw C_t value less than 35 indicates the presence of secondary amplification products. See the Troubleshooting section of this User Manual.
 - c. No Template Control (NTC)
A raw C_t value of 35 or greater indicates a lack of general DNA contamination.
A raw C_t value less than 35 indicates the presence of general DNA contamination in the PCR. See the Troubleshooting section of this User Manual and the Note on Preparing a Workspace Free of DNA Contamination earlier in this protocol.
3. Calculate the ΔC_t for each miRNA in each sample and each dilution.
$$\Delta C_t = C_t^{\text{MOI}} - C_t^{\text{AVG HKG}}$$
4. When biological and/or technical replicates are performed, calculate the average ΔC_t value of each miRNA assay across those replicates for each treatment group.
5. Calculate the $\Delta\Delta C_t$ for each miRNA between your two PCR samples or groups.
$$\Delta\Delta C_t = \Delta C_t (\text{group 2}) - \Delta C_t (\text{group 1})$$

Where group 1 is the control and group 2 is the experimental
6. Calculate the fold-change for each miRNA from group 1 to group 2 as $2^{-\Delta\Delta C_t}$.

OPTIONAL: *If the fold-change is greater than 1, then the result may be reported as a fold up-regulation. If the fold-change is less than 1, then the negative inverse of the result may be reported as a fold down-regulation. The fold-change ratios may also be reported as is.*

NOTE: Detailed Mathematical Explanation of $\Delta\Delta C_t$ Data Analysis Method

Due to the inverse proportional relationship between the threshold cycle (C_t) and the original miRNA expression level, and the doubling of the amount of product with every cycle, the original expression level (L) for each miRNA of interest is expressed as:

$$L = 2^{-C_t}$$

To normalize the expression level of a miRNA of interest (GOI) to a housekeeping sequence (HKG), the expression levels of the two miRNA sequences are divided:

$$\frac{2^{-C_t(\text{GOI})}}{2^{-C_t(\text{HKG})}} = 2^{-[C_t(\text{GOI}) - C_t(\text{HKG})]} = 2^{-\Delta C_t}$$

To determine fold change in miRNA expression, the normalized expression of the GOI in the experimental sample is divided by the normalized expression of the same GOI in the control sample:

$$\frac{2^{-\Delta C_t(\text{expt})}}{2^{-\Delta C_t(\text{control})}} = 2^{-\Delta\Delta C_t} \quad \text{Where } \Delta\Delta C_t \text{ is equal to } \Delta C_t(\text{expt}) - \Delta C_t(\text{control})$$

The complete calculation is as follows:

$$\frac{\frac{2^{-\Delta C_t(\text{GOI}) \text{ expt}}}{2^{-\Delta C_t(\text{HKG}) \text{ expt}}}}{\frac{2^{-\Delta C_t(\text{GOI}) \text{ control}}}{2^{-\Delta C_t(\text{HKG}) \text{ control}}}} = \frac{2^{-[C_t(\text{GOI}) - C_t(\text{HKG})] \text{ expt}}}{2^{-[C_t(\text{GOI}) - C_t(\text{HKG})] \text{ control}}} = \frac{2^{-\Delta C_t \text{ expt}}}{2^{-\Delta C_t \text{ control}}} = 2^{-\Delta\Delta C_t}$$

E. Data Analysis: Optional Standard Curve Method

1. Raw Data

Determine the threshold cycle (C_t) value for each reaction using your instrument's software.

2. Generate a standard curve for each MOI and each HK assay

Plot C_t values of the assays using synthetic miRNA templates on the y-axis versus their dilution factors in a log-scale on the x-axis. Fit the data to a straight line. The x-axis value is directly related to the relative expression level (L) of the miRNA sequence.

3. Determine Relative Expression of Unknowns

Use the standard curve or linear fit equation to convert C_t values to relative expression levels for each MOI or HK sequence in each sample. Be sure that the C_t values fall within the linear range of the appropriate standard curve.

4. Normalization

Divide each MOI relative expression levels by the corresponding HK relative expression levels. Use MOI and HK data from the same dilution of the same sample.

5. Fold-Changes

To determine the fold-change in expression for each miRNA sequence of interest between two different samples, calculate the ratio of its normalized expression levels (again determined from the same PCR conditions) between those samples.

VI. Troubleshooting and FAQs

A. Appearance of multiple PCR products (dissociation curve peaks):

1. Assay Primers Amplify from Non-miRNA cDNA:

If total RNA, instead of small RNA, is used in the miRNA qPCR, more non-specific products and higher background are expected. In this case, using small RNA will normally resolve the problem. Be sure to use the miRNeasy Mini Kit (cat. no. 217004) to enrich for small RNA from your biological samples of total RNA samples, previously purified with the TRIzol[®] procedure.

If the raw C_t value from the assay is greater than 35, consider the miRNA as “Not Detectable” and discounted.

B. Real-time PCR C_t values are too high (> 35 or not detectable):

1. Experimental error:

Use cDNA template generated from the synthetic miRNA sequence of interest as a positive control to troubleshoot the PCR reagents and experimental procedure.

2. Poor RNA quality:

Be sure to perform all recommended quality control checks on the RNA sample. Poor quality RNA can inhibit enzyme activity during reverse transcription generating an insufficient amount of template during the reverse transcription reaction.

3. Not enough template:

- a. Use more input RNA for reverse transcription especially if the lower end of the recommended range had been used previously.
- b. Use a larger volume of template per reaction, but do not use more than 2.0 µl of template per 25-µl reaction. Remember to use the same volume of template in each reaction.

C. Real-time PCR C_t values are too low (< 12): Too much template

1. Use less input RNA for reverse transcription especially if the higher end of the recommended range had been used previously.
2. Use a smaller volume of template per reaction or use diluted template, but do not use less than 1 µl of template per each 25-µl reaction. Remember to use the same volume of template in each reaction.

D. Expression is seen when it is not expected OR The No Reverse Transcription Control (NRT) yields a C_t value < 35:

If total RNA, instead of small RNA, is used in the miRNA qPCR, more non-specific products and higher background are expected. In this case, using small RNA will normally resolve the problem. Be sure to use the miRNeasy Mini Kit (cat. no. 217004) to enrich for small RNA from your biological samples of total RNA samples, previously purified with the TRIzol procedure.

E. The No Template Control (NTC) yields a C_t value < 35:

1. DNA contamination of other reagents, tips, and tubes:
See the Note on Preparing a Workspace Free of DNA Contamination at the beginning of the protocol in this Handbook.
2. Primer Dimers:
Verify presence of primer dimers (< 50 bp in size) by agarose gel electrophoresis. Be sure to use the appropriate RT² SYBR Green qPCR Mastermix to prevent the appearance of primer dimers.

B. Frequently Asked Questions:

1. Will pipetting error affect the miRNA qPCR Assay results?

The passive reference dyes in the PCR master mixes, such as ROX and Fluorescein, are used by the real-time PCR systems to normalize variation from well to well. Therefore, these systems tolerate volume variations caused by pipetting error and evaporation.

2. How can I prevent the evaporation of reaction volume from the wells?

Be sure to carefully and completely seal the PCR plate with the optical thin-wall 8-cap strips or the optical adhesive film before placing it into your thermal cycler. Also, be sure to use a compression pad when using an optical film seal as directed by the manufacturer of your real-time PCR instrument.

3. How reliable are the results from the miRNA qPCR Assays?

Assuming the use of good, consistent experimental technique, real-time PCR methods such as the miRNA qPCR Assays provide very reproducible results. To insure the reliability of your results and to reliably detect smaller fold changes in miRNA expression from the PCR Assays, the performance of replicate determinations (such as biological triplicates) is highly recommended.

If you have additional questions, please check our website (www.SABiosciences.com) for a more complete listing of Frequently Asked Questions (FAQs), or call our Technical Support Representatives at 1-888-503-3187 or 301-682-9200.

Ordering Information

Product	Contents	Cat. no.
RT ² miRNA qPCR Assay (200)	miRNA-specific primer and universal primer for 200 miRNA quantification reactions	Varies

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Notes

RT² miRNA qPCR Assays

Notes

Trademarks: QIAGEN®, QIAzol® (QIAGEN Group); SYBR® (Molecular Probes, Inc.); Roche®, LightCycler® (Roche Group); ROX™ (Applied Biosystems Corporation or its subsidiaries); Eppendorf®, Mastercycler® (Eppendorf AG); Stratagene® (Stratagene); Bio-Rad®, iCycler®, Chromo4™, DNA Engine Opticon®, iQ™5, MyiQ™ (Bio-Rad Laboratories, Inc.); Agilent® (Agilent Technologies, Inc.); LabChip® (Caliper Technologies Corp.); TRIZol® (Molecular Research Center, Inc.).

Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

Limited License Agreement

Use of this product signifies the agreement of any purchaser or user of the RT² miRNA qPCR Assay to the following terms:

1. The RT² miRNA qPCR Assay may be used solely in accordance with the *RT² miRNA qPCR Assay Handbook* and for use with components contained in the Kit only. QIAGEN grants no license under any of its intellectual property to use or incorporate the enclosed components of this Kit with any components not included within this Kit except as described in the *RT² miRNA qPCR Assay Handbook* and additional protocols available at www.qiagen.com.
2. Other than expressly stated licenses, QIAGEN makes no warranty that this Kit and/or its use(s) do not infringe the rights of third-parties.
3. This Kit and its components are licensed for one-time use and may not be reused, refurbished, or resold.
4. QIAGEN specifically disclaims any other licenses, expressed or implied other than those expressly stated.
5. The purchaser and user of the Kit agree not to take or permit anyone else to take any steps that could lead to or facilitate any acts prohibited above. QIAGEN may enforce the prohibitions of this Limited License Agreement in any Court, and shall recover all its investigative and Court costs, including attorney fees, in any action to enforce this Limited License Agreement or any of its intellectual property rights relating to the Kit and/or its components.

For updated license terms, see www.qiagen.com.

PRODUCT WARRANTY

This warranty limits our liability to the replacement of this product in the event the product fails to perform due to any manufacturing defect. SABiosciences Corporation makes no other warranties of any kind, expressed or implied, including without limitation, warranties of merchantability or fitness for a particular purpose. SABiosciences Corporation shall not be liable for any direct, indirect, consequential or incidental damages arising out of the use, the results of use or the inability to use this product.

LIMITED LICENSE STATEMENTS

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 5,079,352, 5,789,224, 5,618,711, 6,127,155, 5,677,152 (claims 1 to 23 only), 5,773,258 (claims 1 and 6 only), 5,407,800, 5,322,770, 5,310,652, 5,994,056, 6,171,785, and claims outside the US corresponding to US Patent No. 4,889,818. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim (such as apparatus or system claims in US Patent No. 6,814,934) and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

This product is provided under an agreement between Molecular Probes, Inc. and SABiosciences and the manufacture, use, sale, or import of this product is subject to one or more of U.S. Patent Nos. 5,436,134; 5,658,751 and corresponding international equivalents, owned by Invitrogen Corp. The purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in research conducted by the buyer, where such research does not include testing, analysis or screening services for any third party in return for compensation on a per test basis. The buyer cannot sell or otherwise transfer (a) this product (b) its components or (c) materials made using this product or its components to a third party or otherwise use this product or its components or materials made using this product or its components for Commercial Purposes. Commercial Purposes means any activity by a party for consideration and may include, but is not limited to: (1) use of the product or its components in manufacturing; (2) use of the product or its components to provide a service, information, or data; (3) use of the product or its components for therapeutic, diagnostic or prophylactic purposes; or (4) resale of the product or its components, whether or not such product or its components are resold for use in research. For information on purchasing a license to this product for purposes other than research, contact Molecular Probes, Inc., Business Development, 29851 Willow Creek Road, Eugene, OR 97402 Tel: (541) 465-8300, Fax: (541) 335-0504.

The purchase of this product includes a limited, non-transferable license under specific claims of U.S. Patent Nos. 6,174,670, 6,569,627 and 5,871,908, owned by the University of Utah Research Foundation or Evotec Biosystems GmbH and licensed to Idaho Technology, Inc. and Roche Diagnostics GmbH, to use only the enclosed amount of product according to the specified protocols. No right is conveyed, expressly, by implication, or by estoppel, to use any instrument or system under any claim of U.S. Patent Nos. 6,174,670, 6,569,627 and 5,871,908, other than for the amount of product contained herein.

Patent Pending.

© 2011 QIAGEN, all rights reserved.

www.qiagen.com

Australia ■ Orders 1-800-243-800 ■ Fax 03-9840-9888 ■ Technical 1-800-243-066

Austria ■ Orders 0800-28-10-10 ■ Fax 0800-28-10-19 ■ Technical 0800-28-10-11

Belgium ■ Orders 0800-79612 ■ Fax 0800-79611 ■ Technical 0800-79556

Brazil ■ Orders 0800-557779 ■ Fax 55-11-5079-4001 ■ Technical 0800-557779

Canada ■ Orders 800-572-9613 ■ Fax 800-713-5951 ■ Technical 800-DNA-PREP (800-362-7737)

China ■ Orders 86-21-3865-3865 ■ Fax 86-21-3865-3965 ■ Technical 800-988-0325

Denmark ■ Orders 80-885945 ■ Fax 80-885944 ■ Technical 80-885942

Finland ■ Orders 0800-914416 ■ Fax 0800-914415 ■ Technical 0800-914413

France ■ Orders 01-60-920-926 ■ Fax 01-60-920-925 ■ Technical 01-60-920-930 ■ Offers 01-60-920-928

Germany ■ Orders 02103-29-12000 ■ Fax 02103-29-22000 ■ Technical 02103-29-12400

Hong Kong ■ Orders 800 933 965 ■ Fax 800 930 439 ■ Technical 800 930 425

Ireland ■ Orders 1800 555 049 ■ Fax 1800 555 048 ■ Technical 1800 555 061

Italy ■ Orders 800-789-544 ■ Fax 02-334304-826 ■ Technical 800-787980

Japan ■ Telephone 03-6890-7300 ■ Fax 03-5547-0818 ■ Technical 03-6890-7300

Korea (South) ■ Orders 080-000-7146 ■ Fax 02-2626-5703 ■ Technical 080-000-7145

Luxembourg ■ Orders 8002-2076 ■ Fax 8002-2073 ■ Technical 8002-2067

Mexico ■ Orders 01-800-7742-639 ■ Fax 01-800-1122-330 ■ Technical 01-800-7742-436

The Netherlands ■ Orders 0800-0229592 ■ Fax 0800-0229593 ■ Technical 0800-0229602

Norway ■ Orders 800-18859 ■ Fax 800-18817 ■ Technical 800-18712

Singapore ■ Orders 1800-742-4362 ■ Fax 65-6854-8184 ■ Technical 1800-742-4368

Spain ■ Orders 91-630-7050 ■ Fax 91-630-5145 ■ Technical 91-630-7050

Sweden ■ Orders 020-790282 ■ Fax 020-790582 ■ Technical 020-798328

Switzerland ■ Orders 055-254-22-11 ■ Fax 055-254-22-13 ■ Technical 055-254-22-12

UK ■ Orders 01293-422-911 ■ Fax 01293-422-922 ■ Technical 01293-422-999

USA ■ Orders 800-426-8157 ■ Fax 800-718-2056 ■ Technical 800-DNA-PREP (800-362-7737)

