

## Release Note: EZ2<sup>®</sup> Connect Protocol Package v8

Dear valued customer,

At QIAGEN<sup>®</sup>, we are continuously making improvements to provide the highest level of performance with our instruments and protocols. Consequently, we are pleased to inform you about the opportunity to upgrade your EZ2 Connect instrument to Protocol Package v8. This protocol package includes the improvements and changes to the protocols listed below.

Protocol name	New protocol version	Changes
EZ1&2 <sup>™</sup> DNA Blood 200 µL Kit	4	<ul style="list-style-type: none"> <li>More accurate run time displayed</li> <li>Implemented additional measures to reduce the risk of lifting the tube up</li> </ul>
EZ1&2 DNA Blood 350 µL Kit	5	<ul style="list-style-type: none"> <li>More accurate run time displayed</li> <li>Sample volume displayed in the loading instructions</li> </ul>
AdnaTest CTC Select Kit	4	<ul style="list-style-type: none"> <li>Implemented additional measures to reduce the risk of heater time out with unstable or high environmental temperature</li> </ul>
EZ1&2 DNA FFPE Kit	3	<ul style="list-style-type: none"> <li>Implemented additional measures to reduce the risk of splashing and mixing issues</li> </ul>
EZ2 RNA FFPE Kit	4	<ul style="list-style-type: none"> <li>Implemented additional measures to reduce the risk of splashing and mixing issues</li> <li>Sample volume displayed in the loading instructions</li> </ul>
EZ2 AllPrep <sup>®</sup> DNA/RNA FFPE Kit	4	<ul style="list-style-type: none"> <li>Implemented additional measures to reduce the risk of splashing and mixing issues</li> <li>More flexible choice of elution volumes. For more information, please refer to "Important Note: EZ2 AllPrep DNA/RNA FFPE Kit and EZ1&amp;2 ccfDNA Kit".</li> </ul>
EZ1&2 ccfDNA Kit	5	<ul style="list-style-type: none"> <li>Implemented additional measures to reduce the risk of heater time out with unstable or high environmental temperature</li> <li>More flexible choice of elution volumes. For more information, please refer to "Important Note: EZ2 AllPrep DNA/RNA FFPE Kit and EZ1&amp;2 ccfDNA Kit".</li> <li>New protocol for urine as starting material. For more information, please refer to "Important Note: EZ2 AllPrep DNA/RNA FFPE Kit and EZ1&amp;2 ccfDNA Kit".</li> </ul>

If you have further questions, please contact your local QIAGEN representative or contact our Technical Support Center at [www.qiagen.com/support/technical-support](http://www.qiagen.com/support/technical-support).

Best regards,

Your EZ2 Connect Team

Trademarks: QIAGEN<sup>®</sup>, Sample to Insight<sup>®</sup>, AllPrep<sup>®</sup>, EZ1&2<sup>™</sup>, EZ2<sup>®</sup> (QIAGEN Group). Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

PROM-22313-002 08/2023 © 2023 QIAGEN, all rights reserved.