

RT² Profiler PCR Arrays

For pathway-focused gene expression analysis

RT² Profiler PCR Arrays are a highly reliable, sensitive gene expression profiling real-time PCR-based technology for analyzing focused panels of genes involved in biological processes, signal transduction, or disease research pathways.

RT² Profiler PCR Arrays provide:

- Focused gene expression profiling for any lab with a real-time instrument
- Integrated controls that assess sample quality and PCR process
- Free easy-to-use data analysis tools (Web- or Excel®-based)

With RT² Profiler PCR Arrays, you can analyze the expression of the most relevant genes in your pathway of interest with complete confidence in your results, to build a better biological story. The RT² Profiler PCR Array system comprises a complete experimental workflow, from sample preparation through data analysis and interpretation, starting with just 25 ng of RNA (or <1 ng of RNA with the RT² PreAMP cDNA Synthesis Kit) and provides free data analysis software. Each catalogued RT² Profiler PCR Array incorporates laboratory-verified assays for 84 pathway-focused genes, 5 housekeeping genes for normalization, and controls that check for sample quality and reaction quality (Figure 1).

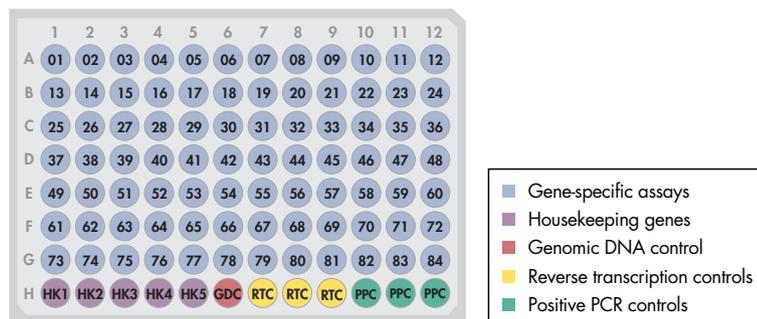


Figure 1. Each well in an RT² Profiler PCR Array measures the relative expression of a gene related to a pathway or disease state. Arrays are available in 4 formats: the 96-well plate, the 384-well plate, the 100-well disc, and the 96x96 chip. Every format includes comprehensive, proprietary controls in addition to the laboratory-verified gene assays, including controls to monitor genomic DNA contamination (GDC), first strand synthesis (RTC), and real-time PCR efficiency (PPC).

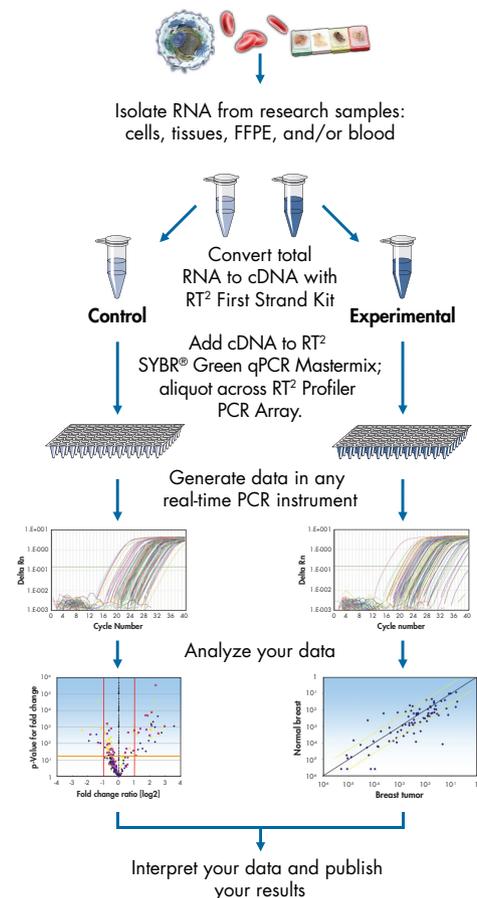


Figure 2. RT² Profiler PCR Array workflow. Starting from cell, tissue, blood, formalin-fixed paraffin-embedded (FFPE), or laser-capture microdissection (LCM) samples, a comprehensive gene expression profile for your pathway of interest can be achieved in less than 3 hours.

Profiling of sample types:

- Cells
- Tissues
- Blood
- FFPE
- LCM



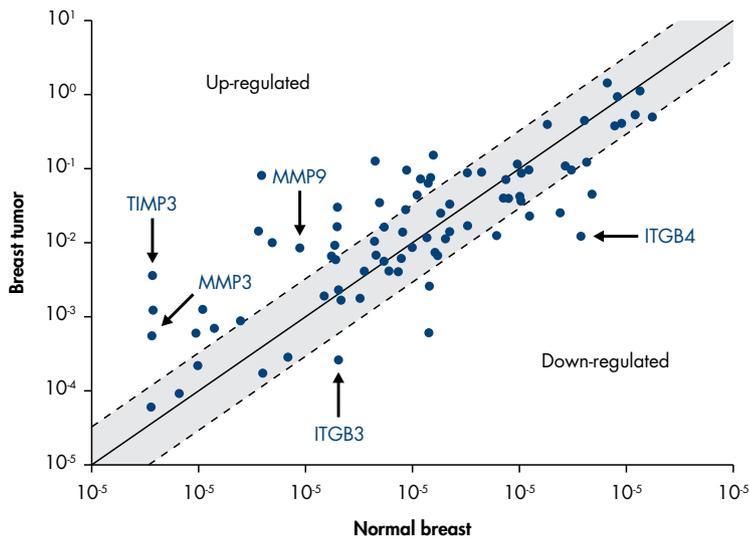


Figure 3. Identification of potential breast cancer biomarkers through expression analysis of extracellular matrix genes. Triplicate total RNA samples from normal human breast and a human breast tumor were characterized using the Human Extracellular Matrix and Adhesion Molecules RT² Profiler PCR Array. Genes encoding the matrix metalloproteinases (MMP3 and MMP9) and their inhibitors (TIMP3) are up-regulated, while genes encoding integrins (ITGB3 and ITGB4) are down-regulated by at least 3-fold in breast tumors relative to normal tissue.

Popular RT² Profiler PCR Arrays for research

Angiogenesis	Inflammatory Cytokines & Receptors
Apoptosis	Innate & Adaptive Immune Response
Autophagy	Mitochondrial Energy Metabolism
Cancer PathwayFinder	NFκB Signaling Pathway
Cell Cycle	Oxidative Stress
Epigenetic Chromatin Modification Enzymes	Signal Transduction PathwayFinder
Epithelial to Mesenchymal Transition (EMT)	TGFβ/BMP Signaling Pathway
Extracellular Matrix & Adhesion Molecules (Figure 3)	Toll-Like Receptor Signaling Pathway
Fibrosis	WNT Signaling Pathway

Over 140 more pathways available online:
www.sabiosciences.com/ArrayList.php

Ordering Information

Product	Contents	Cat. no.
RT ² Profiler PCR Arrays	Pathway, disease, or custom* panels of gene assays	330231
RT ² SYBR Green qPCR Mastermixes	Reagents for real-time PCR reactions (available with ROX, fluorescein, or no internal reference dye)	Varies
RT ² First Strand Kit (12)	Reagents for cDNA synthesis reactions for 12 samples	330401
RNeasy Mini Kit (50)**	Columns, plasticware, and reagents for 50 preps	74104
RNase-free DNase Set (50)	Reagents for 50 preps	79254

Compatible with real-time PCR instruments from the following manufacturers:

QIAGEN®, Agilent/Stratagene®, Applied Biosystems®, Bio-Rad®, Eppendorf®, Fluidigm®, Roche®, TaKaRa®

* For custom arrays, please visit www.sabiosciences.com/custompcrplate.php

** For cell samples. For other sample types, please visit QIAGEN.com

Find out more, visit www.sabiosciences.com/RT2PCR.php !

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor.

Trademarks: QIAGEN® (QIAGEN Group); Applied Biosystems®, SYBR® (Life Technologies); Bio-Rad® (Bio-Rad Laboratories, Inc.); Eppendorf® (Eppendorf AG); Excel® (Microsoft Corporation); Fluidigm® (Fluidigm Corp.); Roche® (Roche Group); Stratagene® (Agilent Technologies Inc.); TaKaRa® (TaKaRa Bio Inc.)
 1071185 1/2012 © 2012 QIAGEN, all rights reserved.

www.SABiosciences.com
www.qiagen.com
 USA = 1-888-503-3187
 Australia = 1-800-243-066
 Austria = 00800-22448000
 Belgium = 00800-22448000

Brazil = 0800-557779
 Canada = 0800-362-7737
 China = 0800-988-0325
 Denmark = 00800-22448000
 Finland = 00800-22448000
 France = 00800-22448000

Germany = 00800-22448000
 Ireland = 00800-22448000
 Italy = 00800-22448000
 Japan = 03-5632-9610
 Luxembourg = 00800-22448000
 Mexico = 01-800-7742-436

Netherlands = 00800-22448000
 Norway = 00800-22448000
 Singapore = 1800-742-4368
 Sweden = 00800-22448000
 Switzerland = 00800-22448000
 UK = 00800-22448000

